

BIOBASIC

INSTRUCTIONS FOR USE

Models: 210, 310, 410 & 600

Before you proceed

This instructions for use is intended for the following product series:

BIOBASIC

We recommend that you read this instructions for use through thoroughly before using the cabinet for the first time. Gram Commercial does not guarantee safe operation if the cabinet is used for anything other than its intended use. Contents of the instructions for use can be subject to change without notice. No part of this instructions for use may be reproduced in any form without expressed written consent of Gram Commercial. Gram Commercial guarantees the cabinet under certain warranty conditions. Gram Commercial is not responsible for any loss or damage of content.

This instructions for use should be considered an integral part of the cabinet and should be stored close to the cabinet and be easy to access. If the instructions for use is lost, please refer to your local distributor or Gram Commercial for a replacement. For current versions of the instructions for use, please go to **www.gram-bioline.com**.

Gram BioLine equipment is designed to be used in a system with monitored additional independent alarms to ensure timely reaction to alarms and thereby maximum item safety



In the event of need for product support, do not hesitate to contact us at:
support@gram-bioline.com

In the event of technical difficulties always contact Gram BioLine technical support or a Gram BioLine authorized service partner.

Copyright © 2006-2022 Gram BioLine, a division of Gram Commercial, Denmark. All rights reserved. The content of this publication is owned by Gram BioLine, unless otherwise noted, and is protected by Danish and international copyright laws and provisions. Information and images may not be used, copied or transferred without the express written permission by Gram BioLine.

Table of Content

Before you proceed	2
Table of Content	4
Intended use	6
Temperature setpoint range and ambient requirements	6
Symbols used	7
Installation	8
Initial setup steps	8
Adjusting the base	9
Anti tilt bracket	10
Wall mounting	11
Surroundings	12
Product inspection	13
Voltage free contact	14
Connecting to power	16
Equipotential bonding	20
Start up	22
The digital display controls	22
General introduction to the controller interface	23
Alarm settings	24
Buttons and useful short cuts	25
Operation parameters	26
Error codes	27
Latching alarms: A2, A3	27
Sensor offset	28
Ordinary use	30
Door gaskets	32
General info	33
Responsibility	33
General info	34
Service	34

Type number plate	35
Access port	36
Regular maintenance	38
Cleaning.....	38
IMPORTANT	40
Disposal.....	41
Datasheet	42
General data: BIOBASIC 210, 310, 410	42
BIOBASIC RR210 – with solid door	44
BIOBASIC RR210 – with glass door	45
BIOBASIC RR310 – with solid door	46
BIOBASIC RR310 – with glass door	47
BIOBASIC RR410 – with solid door	48
BIOBASIC RR410 – with glass door	49
BIOBASIC RF210	50
BIOBASIC RF310	51
BIOBASIC RF410	52
General data: BIOBASIC 600	54
BIOBASIC RR600	56
BIOBASIC RF600	57
Declaration of Conformity	58
Wiring diagrams	59
BIOBASIC RR/RF210, RR/RF310, RR/RF410	59
BIOBASIC RR600	60
BIOBASIC RF600	61
Piping diagram	62
Notes	64



Intended use

BIOBASIC refrigerators (RR) and freezers (RF), are designed and manufactured to provide general purpose storage of non-critical laboratory items.

The BIOBASIC range complies with EN/IEC 60079-15, covering electrical apparatus in Category 3, Zone 2 locations where explosive atmospheres may be present.

Enabling placement of BIOBASIC refrigerators and freezers, in Zone 2 areas categorised according to EN/IEC 60079-14.

The user must ensure that the cabinet is used in accordance with its intended use. Abnormal use or use conflicting with the intended use or guidelines stipulated in the product documentation can lead to; danger to patient safety, damage to stored items, damage to cabinet, danger to user. Gram BIOBASIC equipment is designed to be used in a system with monitored additional independent alarms to ensure timely reaction to alarms and thereby maximum item safety.

Temperature setpoint range and ambient requirements

Model + temperature setpoint range	Minimum ambient operating temperature	Maximum ambient operating temperature
BIOBASIC 210, 310, 410		
RR with solid door: +2/+15°C	+10°C	+35°C
RR with glass door: +2/+15°C	+10°C	+32°C
RF: -25/-5°C	+10°C	+35°C
BIOBASIC 600		
RR: +2/+15°C	+10°C	+43°C
RF: -25/-5°C	+10°C	+43°C

Symbols used



Hazard



Risk of electric shock



Risk of material damage



Risk of personal injury



Risk of burning / freezing



Risk of fire/flammable materials



Risk of explosion/explosive materials



Info



ATEX information

Installation

Initial setup steps




Due to safety and operating considerations, the cabinet must not be used outdoors. The cabinet should be installed in a dry and sufficiently ventilated area. To ensure efficient operation, the cabinet should not be installed in direct sunlight or close to heat sources.



The cabinet interior must not be exposed to corrosive atmospheres.



Avoid placement of the cabinet in a chloric/acidic environment due to risk of corrosion.

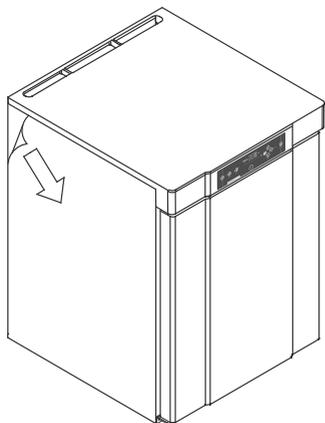

WARNING - POTENTIAL ELECTROSTATIC CHARGING HAZARD

Removing protective packaging and film may cause electrostatic discharge. protective packaging and film shall not be removed in ATEX zones.

The cabinet is shipped with a protective film that shall be removed prior to use.





Clean the cabinet with a mild soap solution prior to use.




The cabinet is only allowed to lay down for very short durations (for instance handling through a doorway). If the cabinet has been laying down, the cabinet must stand up-right for at least 24 hours prior to use. This enables oil in the compressors to run back into place.



Adjusting the base

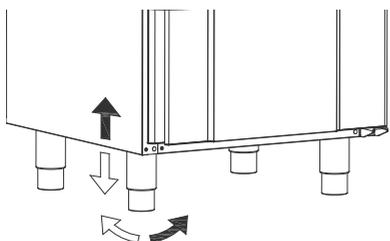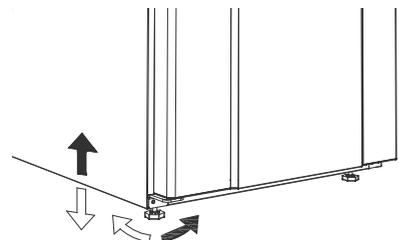

Cabinets equipped with legs should be levelled as shown in the illustrations to the left.

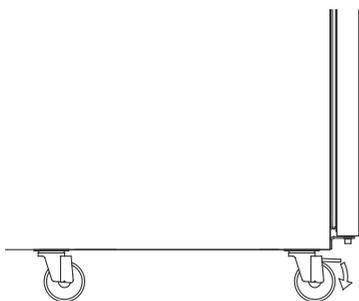

For cabinets equipped with castors, the floor must be level to ensure stable positioning and safe use. When the cabinet is positioned, the 2 front castors should be locked.


Anti tilt bracket



Cabinets with drawers and/or glass door must be secured to a stable vertical surface, ensuring that the cabinet cannot tip over when the drawers are drawn to the outermost position, or the door is open. Brackets for securing the cabinet are included from the factory.



Find the instructions for the tilt bracket below.

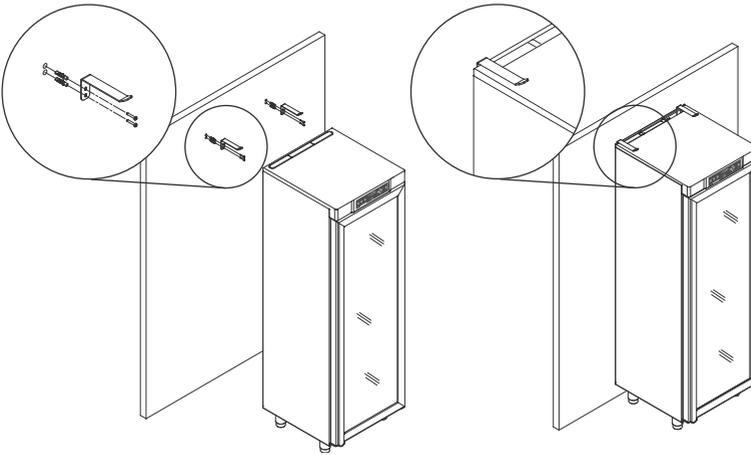

The anti-tilt brackets must be fitted when installing the cabinet, ensuring that the users, surroundings and stored items are not damaged.




Wall mounting

Wall mounting brackets can be supplied if specified. Allowing the cabinet to be lifted from the floor.

Find instructions on wall mounting of a BIOBASIC 210 below, the same procedure applies for mounting 310 and 410.


NB: Applies for 210/310/410 models.

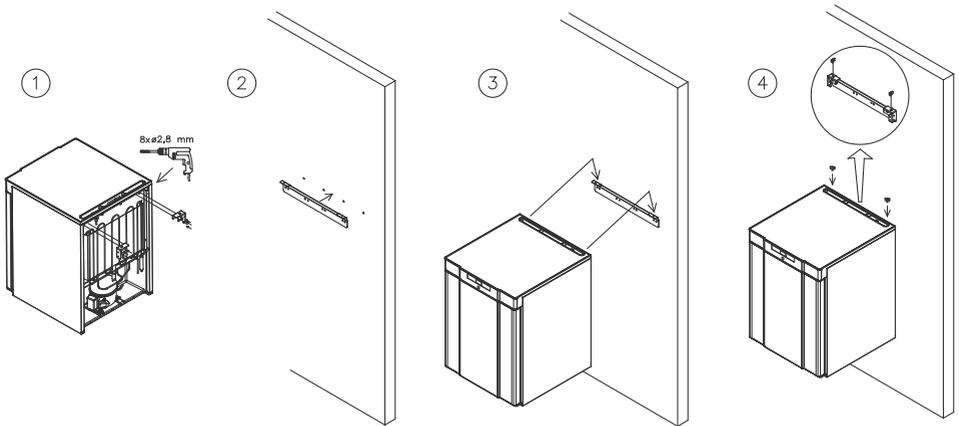

Surroundings



The cabinet must be installed in accordance with the illustrations below.

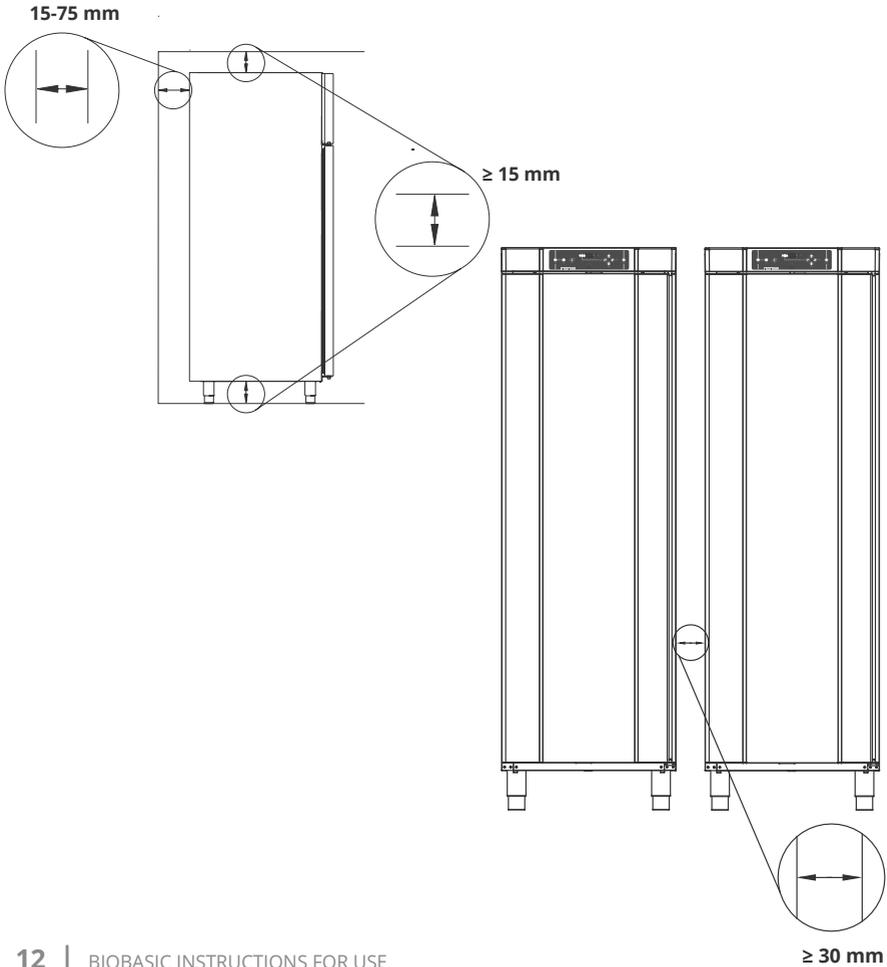



Product inspection



A visual inspection of the cabinet must be conducted prior to putting the cabinet into service.


Check the cabinets structural integrity, that door frames and doors don't have deformities, gaskets seal properly and the doors sit flush up against the door frame.

Voltage free contact


The illustration below shows the three connectors for the relay (e.g. in connecting to CTS or other external monitoring systems). The three connections are respectively Common, NO (Normally Open) and NC (Normally Closed).

The moment voltage is applied, the controller draws the relay, this makes it possible for the controller to respond to both high and low temperature alarms, door alarms and power failures.


Find instructions on setting the alarms in the controller settings section.

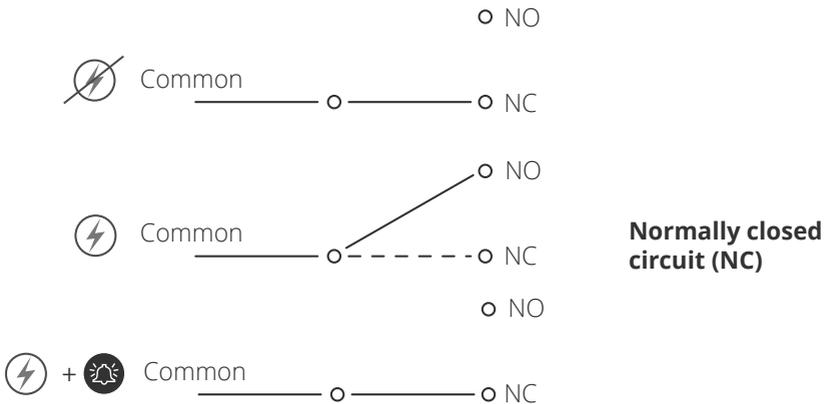

Location of the voltage free contact or 210/310/410

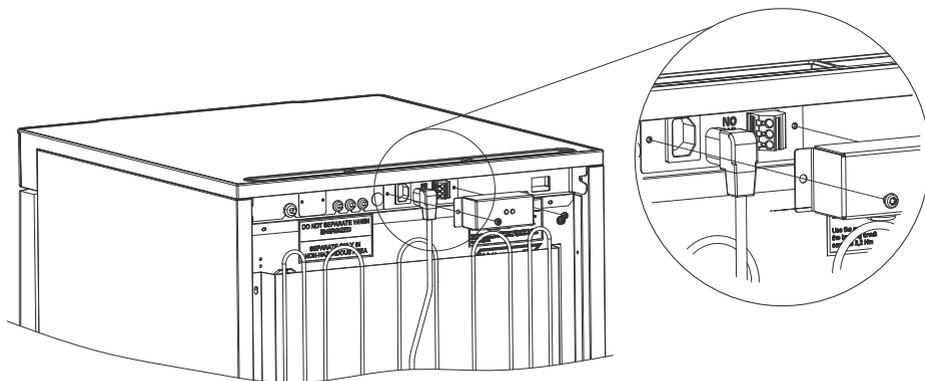

Location of the voltage free contact for 600

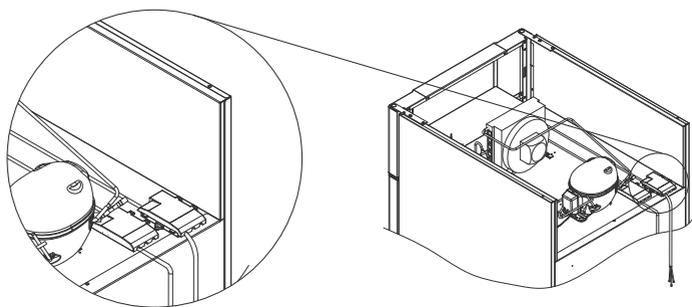

Connection of the voltage free contact should be done by a qualified installer.

Connecting to power



When setting up in an ordinary scenario that is not subject to regulations for EN 60079-15 zone 2:

The appliance may be connected in accordance with applicable local heavy current regulations.


Note that there are special regulations for products that are in accordance with EN 60079-15 zone 2 and EN 60079-14: Explosive atmospheres – Electrical installations design, selection and erection.

The appliance has been manufactured in accordance with EN 60079-15: Electrical apparatus for explosive gas atmospheres – Part 15: Type of protection II 3G Ex ec nC ic IIB Tx Gc. Zone 2 is the applicable zone. If the appliance is to be installed in a zone 2 environment, specialist personnel should perform the installation, or be consulted beforehand, in order to ensure that the appliance is installed in compliance with the guidelines currently contained in the standard. The cabinet is intended for connection to alternating current. The connection values for voltage (V) and frequency (Hz) are given on the type/number plate.




BIOBASIC 210, 310, 410 – The mains terminal is accessed via the preload cover on the back of the cabinet. Unscrew the preload cover to access the mains supply terminal. Ensure that the preload cover is reinstalled after plugging in the power cord. The leaf spring in the preload cover must engage and preload the plug of the cord as shown in the illustrations below.

BIOBASIC 600 – The power cord is plugged in the terminal box on the back of the cabinet. The plug is then fixated in place by the hanger that is built into the terminal box. Please note that the hanger should be fitted tightly around the plug, as shown. In all cases, ensure that the mains plug is seated completely in the terminal of terminal on the cabinet.

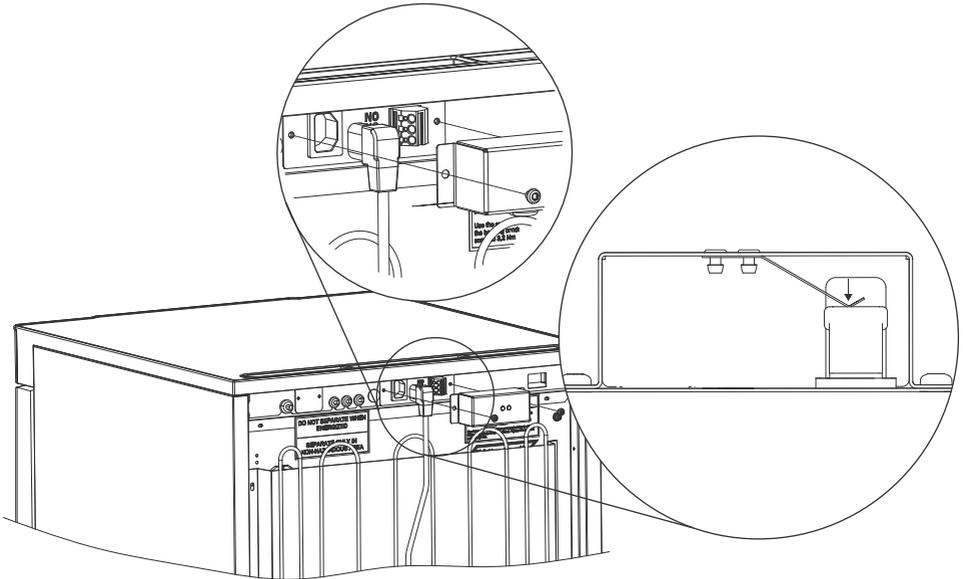

Power cord plugged into 210/310/410 cabinets.

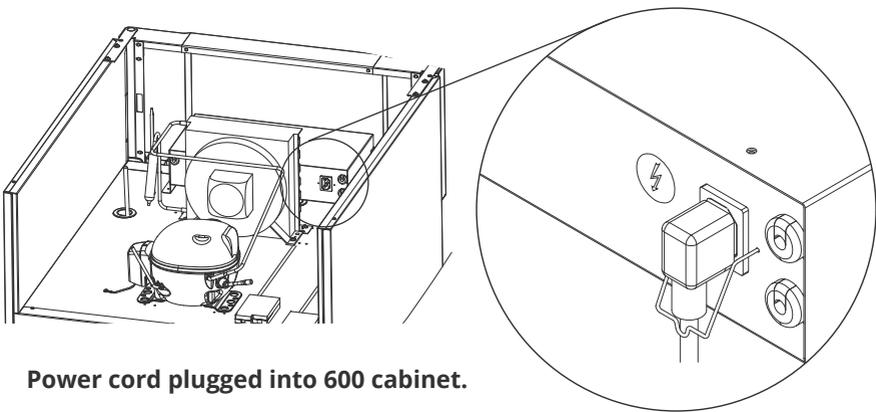

Power cord plugged into 600 cabinet.

The appliance must be connected to the external power supply using a suitable device which mechanically prevents the plug and socket from being separated unintentionally. The connection must be labelled:

“DO NOT SEPARATE WHEN ENERGIZED”.

Fuses and similar must never be removed or replaced while the appliance is connected to a power source. The electrical terminal box must never be opened while the appliance is connected to a power source. The compressor starting equipment must never be dismantled while the appliance is connected to a power source. The LED lighting must never be dismantled while the appliance is connected to a power source. Whenever electrical components are dismantled or replaced, the appliance must be moved to an area in which there is no risk of ignition caused by the electrical components or gases contained in the appliance. Never use the cabinet if the plug is damaged. The cabinet should be examined by a Gram BioLine authorised service technician in such cases.

Please refer to EN 60079-14: Explosive atmospheres – Electrical installations design, selection and erection for installation requirements in an ATEX environment.

In both cases:

Use a three-wire plug, if the power outlet is intended for a three-wire plug, the lead in green/yellow insulation should be connected to the ground terminal. Power must be connected via a wall socket. The wall socket should be easily accessible. All earthing requirements stipulated by the local electricity authorities must be observed. The cabinet plug and wall socket should then give correct earthing. If in doubt, contact your local supplier or authorized electrician.

**Technical support -**

In the event of technical difficulties always contact Gram BioLine technical support or a Gram BioLine authorized service partner.

**For Ex environments -**

Special conditions for safe use may apply to this product when installing in an EN 60079-14 environment.
Please see corresponding Ex certificate for specifications.

**For Ex environments -**

Open containers inside the storage chamber may impact the ATEX zone classification.

Equipotential bonding



For installation in ATEX Cat. 3 Zone 2 areas, it is mandatory to have a equipotential bonding, it is not sufficient to use protective earth through the mains connection.

To secure equipotential bonding of the unit – the mounted external bonding conductor must be used in accordance with national installation requirements e.g. EN 60079-14.


- Mounting of the bonding conductor should be done according to the following illustrations.
- Please find location for connection facilities on the back of the cabinet marked with “Attention – Equipotential bonding”.
- The bonding conductor should be at least 4 mm² gauge.
- Use a ring terminal to ensure adequate bonding.
- Use the supplied M5 machine screw and washer to attach the bonding conductor to the cabinet. Tighten the machine screw to 3.2 Nm.

Bonding of the cabinet is illustrated on the following page.




ATTENTION -

Please note, this location is the only manufacturer approved location for Equipotential bonding.

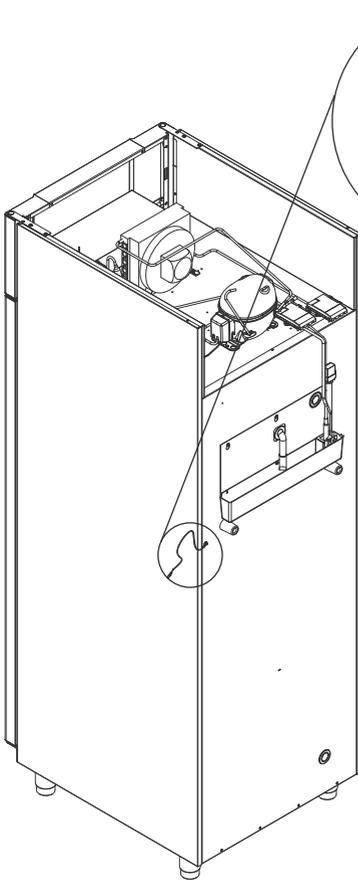

BIOBASIC 600


3.2 Nm

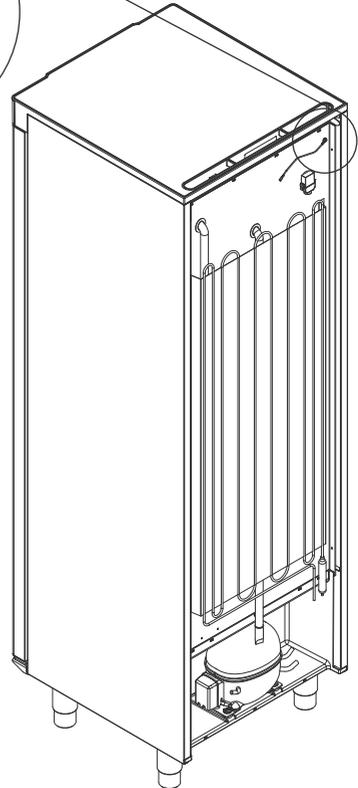

BIOBASIC 210/310/410

Start up

The digital display controls

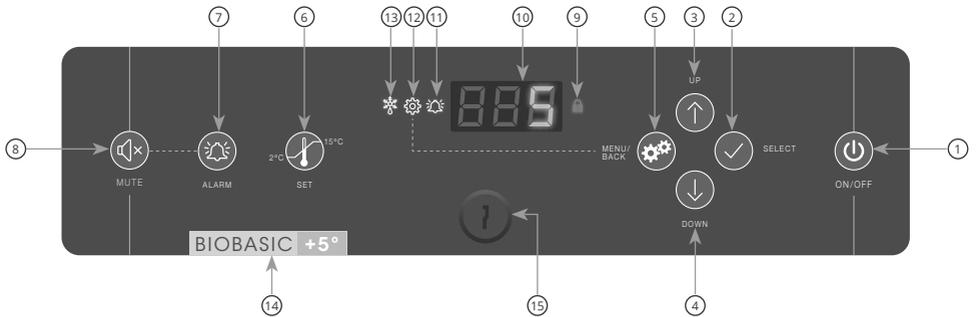

- ① On/Off
- ② Select or confirm a menu parameter
- ③ Navigate upwards in a given menu/raise a given value
- ④ Navigate downwards in a given menu/lower a given value
- ⑤ Parameter settings menu/ go a menu-step back
- ⑥ Setpoint temperature setting
- ⑦ Temperature alarm settings
- ⑧ Acknowledge alarm, mute for 5 minutes
- ⑨ Keypad lock engaged
- ⑩ Display
- ⑪ Alarm registered
- ⑫ Parameter settings menu is open
- ⑬ Defrost in progress
- ⑭ Visual distinction between refrigerator or freezer
- ⑮ Door lock



General introduction to the controller interface

Turn On/Off:

Shortly press  to turn the cabinet On and press for 6 seconds to turn the cabinet Off.

Initiation procedure:

Readout of the software version and variant will occur shortly after turning the cabinet on. The cabinet will then automatically start a defrost cycle, and terminate it again after a system check.

The cabinet is ready when the temperature is displayed.

The cabinet will always commence operation when initially connected to a power supply. For instance after a power outage or when plugging the cabinet in for the first time.

Setpoint temperature adjustments:

Temperature adjustments are done by pressing  for 6 seconds, prompting the set point temperature. Adjust the set point temperature by either pressing  or . Confirm the settings by pressing .

Alarm settings

Press  for 6 seconds to enter the alarm settings

Alarm settings 		Unit	Explanation
	HL	[°C]	High temperature alarm limit. Code for activated alarm [A2]
	LL	[°C]	Low temperature alarm limit. Code for activated alarm [A3]
	Hd	[Min.]	Delay of high temperature alarm
	Ld	[Min.]	Delay of low temperature alarm
	dA	On/ Off	Door alarm. Code for activated alarm [A1]. [1=on / 0=Off]
	dAd	[Min.]	Delay of door alarm
	bU	On/ Off	Acoustic signal for alarm codes [A1], [A2] and [A3]. [1=on / 0=Off]



Buttons and useful short cuts

Buttons:	Press for:	Function:
	-	Acknowledge alarm, mute for 5 minutes
	> 6 seconds	Access the alarm settings
	> 6 seconds	Access the parameter settings
	> 6 seconds	Adjust/show setpoint temperature value
	> 3 seconds	Manually start or stop a defrost
	> 6 seconds	Activating/deactivating the keypad lock
	-	Show highest registered temperature spike (since the last reset of alarm and temperature history)
	-	Show lowest registered temperature spike (since the last reset of alarm and temperature history)
	> 3 seconds	Clear and reset alarm and temperature history
	> 6 seconds	Restores factory settings

Operation parameters



ATTENTION -

Changing operation parameters without the expressed consent from Gram BioLine can have unintended implications on performance and potentially void warranty.

Press  for 6 seconds to enter operation parameters.

Operation parameters 	↴	Unit	
	CA	[K]	Offset of A-sensor. Reference sensor for refrigeration and alarm system
	d1		Number of defrosts per 24 hours (4 is factory setting)
	d2	[°C]	Termination temperature in the evaporator during a defrost
	Li	On/Off	*Only for glass door models* - Turn light on or Off
	tEr	 ↴	Relay test/Component test
	tC	[°C]	Test compressor relay
	tF	[Min.]	Test evaporator fan
	td	[Min.]	Test defrost element relay (RF models)
	tL	On/Off	Test
	tA	[Min.]	Test alarm relay (will trip voltage free)
	tdP	On/Off	Test display
Sensor read-out	P-A	[°C]	Test output of A-sensor
	P-B	[°C]	Test output of B-sensor



Error codes

Display code	Explanation
- 0 -	Door is open
A1	Door alarm “dAd” has been activated
A2	High temperature alarm “HL” is or has been activated
A3	Low temperature alarm “LL” is or has been activated
F1	Error on the main cabinet sensor. The refrigeration system will use an emergency program to make the cabinet run. Temperature stability will be affected. Service is required
F2	Error on the evaporator sensor. Service is required



ATTENTION -

High and low temperature alarms set up in the cabinets controller must be accompanied by additional redundant independent external alarms to ensure maximum item safety.

Please, find instructions for connecting the “voltage-free contact” in the “Installation” section.

Latching alarms: A2, A3

Due to the potential implications of alarms, the red alarm bell light will continue to remain on to warn the user of an alarm having occurred. Regardless of acknowledging, the light will remain on until an alarm history reset has been performed (holding  +  for more than 3 seconds), note that this will also reset temperature and alarm history.

Sensor offset

Sensor offset is used in cases where there are deviations in the cabinets actual operation compared to control measurements by independent temperature measuring systems.

The A-sensor is used to manage the cabinets refrigeration system, and is also the reference sensor for the display and alarms.

The A-sensor is offset if the actual temperature in the cabinet does not match the setpoint, despite taking the hysteresis into consideration. Offset of A sensor is named "CA".

Offsetting the A-sensor

- ↳ Press and hold  for more than 6 seconds
- ↳ Press  to select "CA"
- ↳ Press  +  to offset the A-sensor
- ↳ Press  to confirm the set value

The A-sensor is now offset, proceed to other parameters by pressing  navigate by using  or .

- ↳ Leave the user menu by pressing  several times until the cabinet temperature is shown in the display.



Practical example of offsetting:

Example 1 – The temperature in the cabinet is operating colder than the actual setpoint.

With a setpoint of +4°C, the actual temperature inside the cabinet is between +2°C and +4°C. The desired temperature range is between +3°C and +5°C.

This means that “CA”, in this case, should be -1.0 K, so that the refrigeration system stops 1.0 K before and starts 1.0 K later than the setpoint normally otherwise would dictate.

Example 2 – The temperature in the cabinet is operating warmer than the actual setpoint.

With a setpoint of +4°C, the actual temperature inside the cabinet is between +4°C and +6°C. The desired temperature range is between +3°C and +5°C.

This means that “CA”, in this case, should be 1.0 K, so that the refrigeration system stops 1.0 K later and starts 1.0 K earlier than the setpoint normally otherwise would dictate.

Ordinary use




The cabinet is not suited for storing items that emit vapours, as they might corrode the cabinet and its components.



All items in the cabinet that are not encapsulated, or wrapped, should be covered to reduce the risk of corrosion of the cabinet and its components.

The cabinet interior must not be exposed to corrosive atmospheres.

Items should be evenly distributed in the cabinet, with minimum layer thickness/maximum surface. And at the same time, the air should be able to circulate freely between the items.

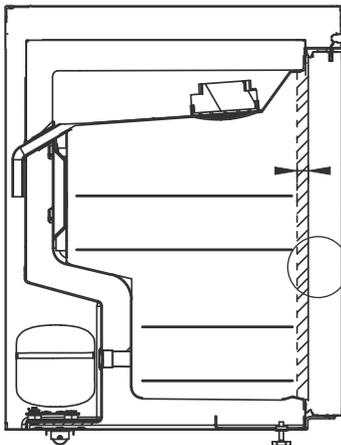


There must be a 20 mm gap between the inner part of the door and the stored items, to enable airflow at the front of the cabinet.




Items placed on the top, bottom or very front of the cabinet will cause the air circulation to be impeded, reducing the cabinet's performance.

Keep the marked areas in the cabinet (shown on this page) clear of all items, ensuring adequate air circulation, and therein cooling.

Do not place items beneath the lowest shelf bracket.

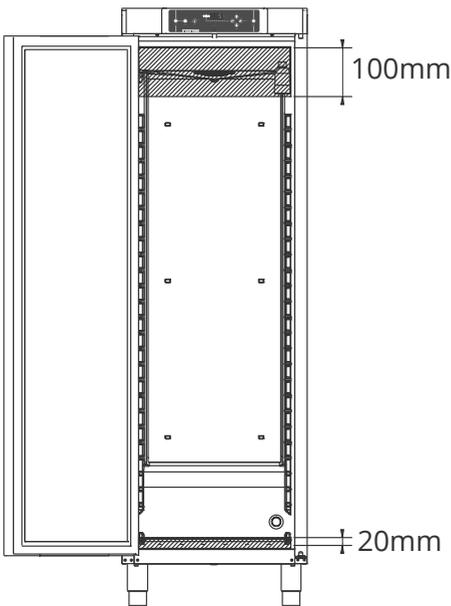

BIOBASIC 210/310/410

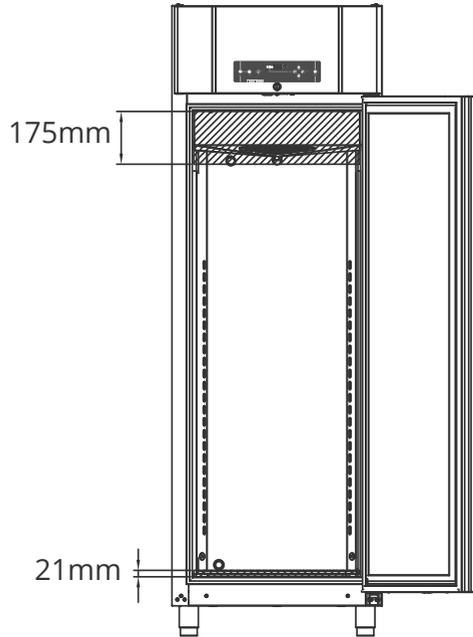


BIOBASIC 600

Door gaskets

Door gaskets are an important part of a cabinet, door gaskets with impaired functionality reduces a cabinets seal with the door.

Impaired seals can lead to increased humidity in the storage chamber, iced evaporator (and thus reduced cooling capacity), and in some cases, decreased lifetime expectancy of the cabinet. It is therefore very important to be aware of the door gaskets condition. Regular inspections are recommended.

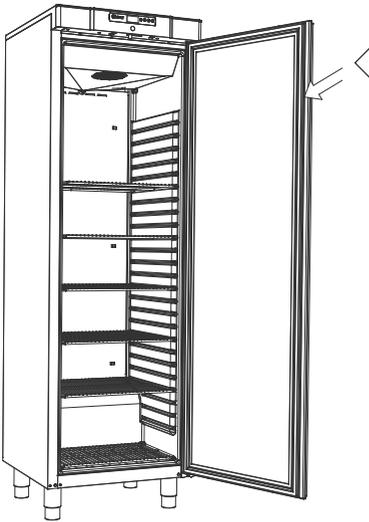

The door gasket should be cleaned regularly with a mild soap solution.

If a gasket is to be replaced, please contact your local Gram BioLine distributor.

General info



Responsibility



Read the following carefully, for information on technical safety and responsibility on Gram BioLine products.

Warranty may be void in the instance the cabinet is being used for applications other than its intended use, or otherwise not in accordance with the guidelines stipulated in the instructions for use.



Defective parts must be replaced with original parts from Gram BioLine. Gram BioLine can only guarantee functional and safety requirements on the cabinets, if above-mentioned is adhered to.



The cabinet should be checked at least once yearly by a Gram BioLine-authorized technician.

General info

Service



WARNING -

DO NOT OPEN, MAINTAIN OR SERVICE IN AN AREA WHEN AN EXPLOSIVE ATMOSPHERE IS PRESENT

The refrigeration system and the hermetically sealed compressor require no maintenance.

However, the condenser and air filter require regular cleaning. If refrigeration fails, first look to see whether the cabinet has been unintentionally switched Off, or whether a fuse has blown.

If the cause of failure cannot be found, contact your supplier quoting Type and S/N, find this on the type number plate.



When servicing -


Make sure the appliance is switched Off at the socket before service is performed on the cabinet. It is not sufficient to switch Off the cabinet on the On/Off  button, as current will persist in some electrical parts of the cabinet.




Be aware that cabinets using hydrocarbons (HC) as refrigerant, may require special handling by qualified technicians.

Type number plate

The type/number-plate located on the inside of the storage chamber, please find reference for location below.

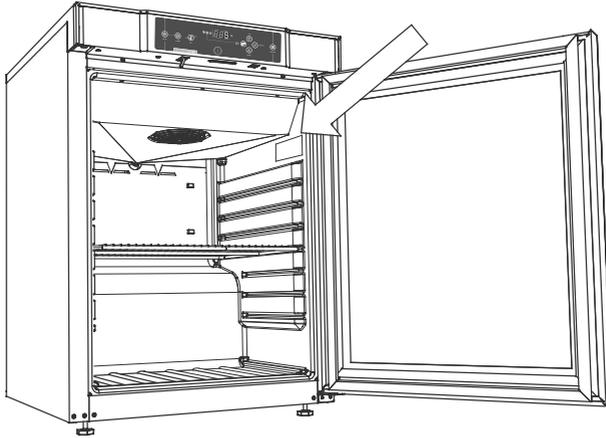

BIOBASIC 210/310/410

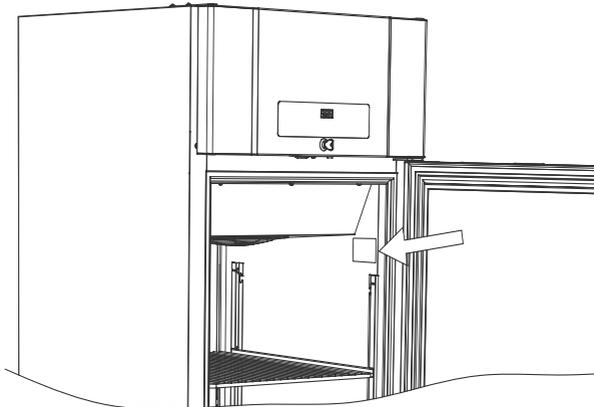


BIOBASIC 600

Access port


All access ports are constructed in the same fashion, with a conical polystyrene plug (1), fitted from the back of the cabinet and a plastic cap (2), fitted from the inside of the cabinet.

Please note, it is very important to re-fit the polystyrene plug and plastic cap after mounting sensor, probe etc. Failing to do so can result in lowered performance or malfunction of the cabinet.

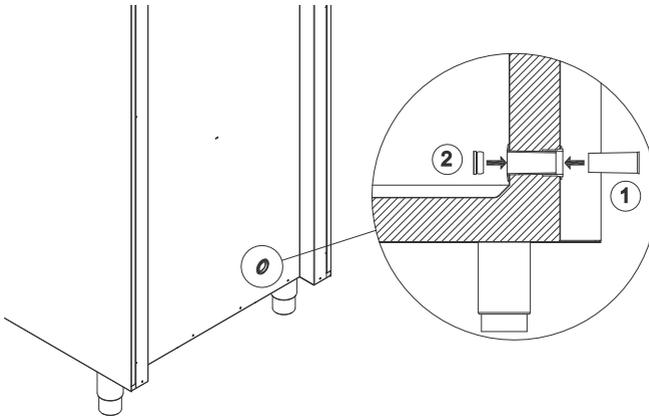




Access ports are clearly marked "Access port" on the cabinet.
Find locations below.

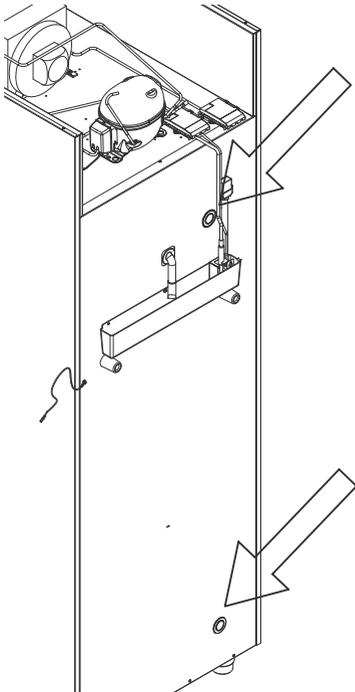

BIOBASIC 600

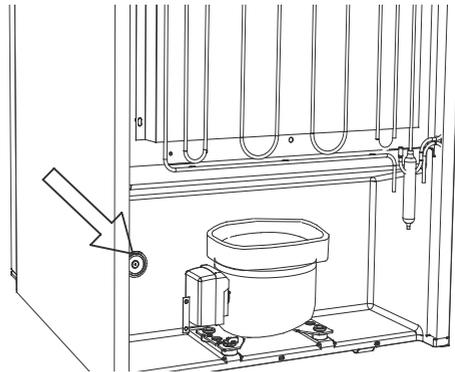


BIOBASIC 210/310/410


Regular maintenance

Cleaning

The cabinet should be cleaned internally with a mild soap solution. (Max. 85°C) at suitable intervals and checked thoroughly before it is put into operation again.

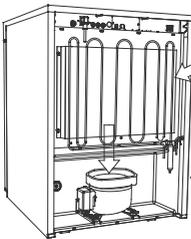

The compressor compartment and in particular the condenser must be kept free from dust and dirt. This is best done with a vacuum cleaner and a brush. Do not flush compressor compartment or evaporator with water as this may cause short circuits in the electrical system.

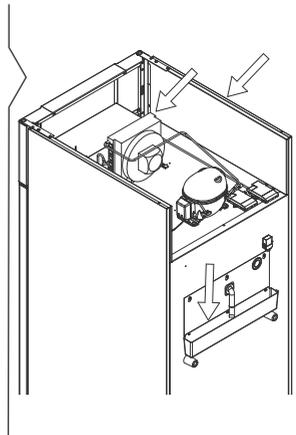

It is recommended that the re-evaporation tray is checked regularly for foreign objects and cleaned accordingly.

Be careful not to damage the defrost water tube and heating element (located in the tray) when cleaning.



Cleaning agents containing chlorine or compounds of chlorine as well as other corrosive agents may not be used, as they might cause corrosion to the stainless panels of the cabinet and the evaporator system.



The cabinet must not be connected to a power source while cleaning.



IMPORTANT



ATTENTION



There may occur sharp edges on the cabinet housing, compressor room, and interior furnishings. Show due diligence when handling the cabinet, neglect of these precautions can lead to injuries.

Danger of wedging in the frame slot between the door and the cabinet, show due diligence when opening and closing the cabinet door. Negligence of these precautions can lead to injuries.

Danger of wedging in the drawer column between the drawers and interior of the cabinet, show due diligence when using the drawers. Negligence of these precautions can lead to injuries.

Unlocked castors can lead to unexpected movements of the cabinet. Lock the castors after installation. Negligence of these precautions can lead to injuries.

The re-evaporation tray, re-evaporation tray heating element, pressure pipes and compressors develop considerable heat during operation. Assure that these components are sufficiently tempered before touching. Negligence of these precautions can lead to injuries.

The evaporator develops considerable cold during operation. Assure that the evaporator is sufficiently tempered before touching. Negligence of this precaution may lead to injuries.

The fan may cause injury during operation, avoid touching the fans while the cabinet is connected to the main power. Negligence of these precautions can lead to injuries.

Do not use electrical appliances inside the cabinet.

Unauthorized modifications to the cabinet can void warranty.



Disposal



Electrical and electronic equipment (EEE) contains materials, components and substances that can be dangerous and harmful to human health and the environment if the waste (WEEE) is not disposed of properly.

Products that are labelled with a “crossed-out wheeled bin” is electric and electronic equipment. The crossed out wheeled bin symbolizes that waste of this type cannot be disposed of with unsorted municipal waste but must be collected separately.

Contact your local Gram BioLine distributor when the cabinet needs to be disposed of.

For additional information, see our website:
www.gram-biobasic.com

Datasheet

General data: BIOBASIC 210, 310, 410

Technical specifications	Data
Connection	230 VAC, 50 Hz
Control Unit	Gram BIOBASIC MPC-46
Alarms	Acoustic and visual temperature and door alarms
Alarm ports	Voltage free contact (230 VAC/ 8A)
Access port	1 pcs. ø24 mm
Door	Right or left hinged
Material interior	PS lining
Materiale exterior	White lacquered steel
Wall thickness	52 mm
Insulation	Polyurethane foam with HFC-free cyclopentane propellant
Air system	Gram BioLine ventilated air distribution system
IP class	IP21



BIOBASIC RR210 – with solid door

Technical specifications	Data
Temperature range	+2°C/+15°C
Ambient temperature range	+10/+35°C
SW variant	K70
K-Value	0.36 W/(m ² *K)
ATEX certificate	DTI 21ATEX0161X
ATEX Marking	II 3G Ex ec nC ic IIB T6 Gc
Gross volume	125 litres
Net volume	104 litres
Dimensions – H x B x D	830 x 595 x 640 mm
Defrost system	Automatic smart defrost with re-evaporation of the defrost water. Air defrost
Refrigerant	R600a
Refrigerant charge	33 g
Refrigeration capacity at -10°C	154 Watt
GWP – CO ₂ e	N/A
Energy consumption – default setpoint	0.61 kWh/24h
Heat emission 100%	97 Watt
Heat emission default setpoint	26 Watt
Nominal consumption	95.1 Watt
Start current	7.6 A
Sound level	36.5 dB(A)

BIOBASIC RR210 – with glass door

Technical specifications	Data
Temperature range	+2°C/+15°C
Ambient temperature range	+10/+32°C
SW variant	K73
K-Value	0.36 W/(m ² *K)
ATEX certificate	DTI 21ATEX0161X
ATEX Marking	II 3G Ex ec nC ic IIB T6 Gc
Gross volume	125 litres
Net volume	104 litres
Dimensions – H x B x D	830 x 595 x 640 mm
Defrost system	Automatic smart defrost with re-evaporation of the defrost water. Air defrost
Refrigerant	R600a
Refrigerant charge	33 g
Refrigeration capacity at -10°C	154 Watt
GWP – CO ₂ e	N/A
Energy consumption – default setpoint	0.61 kWh/24h
Heat emission 100%	102 Watt
Heat emission default setpoint	41 Watt
Nominal consumption	100.1 Watt
Start current	7.6 A
Sound level	36.5 dB(A)

BIOBASIC RR310 – with solid door

Technical specifications	Data
Temperature range	+2°C/+15°C
Ambient temperature range	+10°C/+35°C
SW variant	K70
K-Value	0.36 W/(m ² *K)
ATEX certificate	DTI 21ATEX0161X
ATEX Marking	II 3G Ex ec nC ic IIB T6 Gc
Gross volume	218 litres
Net volume	189 litres
Dimensions – H x B x D	1220 x 595 x 640 mm
Defrost system	Automatic smart defrost with re-evaporation of the defrost water. Air defrost
Refrigerant	R600a
Refrigerant charge	45 g
Refrigeration capacity at -10°C	154 Watt
GWP – CO ₂ e	–
Energy consumption – default setpoint	0.63 kWh/24h
Heat emission 100%	93.7 Watt
Heat emission default setpoint	27.6 Watt
Nominal consumption	99.1 Watt/0.9 A
Start current	7.6 A
Sound level	38.6 dB(A)



BIOBASIC RR310 – with glass door

Technical specifications	Data
Temperature range	+2°C/+15°C
Ambient temperature range	+10°C/+32°C
SW variant	K73
K-Value	0.36 W/(m ² *K)
ATEX certificate	DTI 21ATEX0161X
ATEX Marking	II 3G Ex ec nC ic IIB T6 Gc
Gross volume	218 litres
Net volume	189 litres
Dimensions – H x B x D	1220 x 595 x 640 mm
Defrost system	Automatic smart defrost with re-evaporation of the defrost water. Air defrost
Refrigerant	R600a
Refrigerant charge	45 g
Refrigeration capacity at -10°C	154 Watt
GWP – CO ₂ e	N/A
Energy consumption – default setpoint	0.63 kWh/24h
Heat emission 100%	93.7 Watt
Heat emission default setpoint	27.6 Watt
Nominal consumption	99.1 Watt/0.9 A
Start current	7.6 A
Sound level	38.6 dB(A)

BIOBASIC RR410 – with solid door

Technical specifications	Data
Temperature range	+2°C/+15°C
Ambient temperature range	Solid door +10°C/+35°C
SW variant	K70
K-Value	0.35 W/(m ² *K)
ATEX certificate	DTI 21ATEX0161X
ATEX Marking	II 3G Ex ec nC ic IIB T6 Gc
Gross volume	346 litres
Net volume	312 litres
Dimensions – H x B x D	1876 x 595 x 640 mm
Defrost system	Automatic smart defrost with re-evaporation of the defrost water. Air defrost
Refrigerant	R600a
Refrigerant charge	51 g
Refrigeration capacity at -10°C	207 Watt
GWP – CO ₂ e	N/A
Energy consumption – default setpoint	0.73 kWh/24h
Heat emission 100%	109.6 Watt
Heat emission default setpoint	30.2 Watt
Nominal consumption	113 Watt/0.8 A
Start current	6.4 A
Sound level	35.1 dB(A)



BIOBASIC RR410 – with glass door

Technical specifications	Data
Temperature range	+2°C/+15°C
Ambient temperature range	Glass door +10°C/+32°C
SW variant	K73
K-Value	0.35 W/(m ² *K)
ATEX certificate	DTI 21ATEX0161X
ATEX Marking	II 3G Ex ec nC ic IIB T6 Gc
Gross volume	346 litres
Net volume	312 litres
Dimensions – H x B x D	1876 x 595 x 640 mm
Defrost system	Automatic smart defrost with re-evaporation of the defrost water. Air defrost
Refrigerant	R600a
Refrigerant charge	51 g
Refrigeration capacity at -10°C	207 Watt
GWP – CO ₂ e	N/A
Energy consumption – default setpoint	0.73 kWh/24h
Heat emission 100%	109.6 Watt
Heat emission default setpoint	30.2 Watt
Nominal consumption	113 W/0.8 A
Start current	6.4 A
Sound level	35.1 dB(A)

BIOBASIC RF210

Technical specifications	Data
Temperature range	-25°C/-5°C
Ambient temperature range	+10°C/+35°C
SW variant	F70
K-Value	0.36 W/(m ² *K)
ATEX certificate	DTI 21ATEX0161X
ATEX Marking	II 3G Ex ec nC ic IIB T3 Gc
Gross volume	125 litres
Net volume	104 litres
Dimensions – H x B x D	830 x 595 x 640 mm
Defrost system	Automatic smart defrost with re-evaporation of the defrost water. Heating element
Refrigerant	R600a
Refrigerant charge	30 g
Refrigeration capacity at -25°C	156 Watt
GWP – CO ₂ e	–
Energy consumption – default setpoint	1.1 kWh/24h
Heat emission 100%	145.9 Watt
Heat emission default setpoint	63.9 Watt
Nominal consumption	133 Watt/1.1 A
Start current	6.6 A
Sound level	39.7 dB(A)



BIOBASIC RF310

Technical specifications	Data
Temperature range	-25°C/-5°C
Ambient temperature range	+10°C/+35°C
SW variant	F70
K-Value	0.36 W/(m ² *K)
ATEX certificate	DTI 21ATEX0161X
ATEX Marking	II 3G Ex ec nC ic IIB T3 Gc
Gross volume	218 litres
Net volume	189 litres
Dimensions – H x B x D	1220 x 595 x 640 mm
Defrost system	Automatic smart defrost with re-evaporation of the defrost water. Heating element
Refrigerant	R600a
Refrigerant charge	35 g
Refrigeration capacity at -25°C	204 Watt
GWP – CO ₂ e	N/A
Energy consumption – default setpoint	1.3 kWh/24h
Heat emission 100%	141.5 Watt
Heat emission default setpoint	70.5 Watt
Nominal consumption	152 Watt/1.0 A
Start current	8.0 A
Sound level	39.9 dB(A)

BIOBASIC RF410

Technical specifications	Data
Temperature range	+2°C/+15°C
Ambient temperature range	+10°C/+35°C
SW variant	F70
K-Value	0.35 W/(m ² *K)
ATEX certificate	DTI 21ATEX0161X
ATEX Marking	II 3G Ex ec nC ic IIB T3 Gc
Gross volume	346 litres
Net volume	312 litres
Dimensions – H x B x D	1875 x 595 x 640 mm
Defrost system	Automatic smart defrost with re-evaporation of the defrost water. Heating element
Refrigerant	R600a
Refrigerant charge	47 g
Refrigeration capacity at -25°C	204 Watt
GWP – CO ₂ e	N/A
Energy consumption – default setpoint	1.6 kWh/24h
Heat emission 100%	153.4 Watt
Heat emission default setpoint	88.7 Watt
Nominal consumption	157 Watt/1.0 A
Start current	8.0 A
Sound level	38.6 dB(A)



General data: BIOBASIC 600

Technical specifications	Data
Ambient temperature range	+10°C/+43°C
Connection	230 VAC, 50 Hz
Control Unit	Gram BIOBASIC MPC-46
Alarms	Acoustic and visual temperature and door alarms
Alarm ports	Voltage free contact (230 VAC/8A)
Access port	2 pcs. ø24 mm
Gross volume	610 litres
Net volume	536 litres
Door	Right or left hinged
Material interior	Stainless steel
Materiale exterior	White lacquered steel
Wall thickness	70 mm
K-Value	0.27 W/(m ² *K)
Dimensions – H x B x D	2125 x 700 x 895 mm
Gross weight	45.5 kg
Net weight	42.5 kg
Insulation	Polyurethane foam
Air system	Gram BioLine ventilated air distribution system
IP class	IP21



BIOBASIC RR600

Technical specifications	Data
Temperature range	+2°C/+15°C
SW variant	K70
Defrost system	Automatic smart defrost with re-evaporation of the defrost water. Air defrost
ATEX certificate	DTI 19ATEX0114X
ATEX Marking	II 3G Ex ec nC ic IIB T5 Gc
Refrigerant	R290
Refrigerant charge	75 g
Refrigeration capacity at -10°C	336 Watt
GWP - CO2e	N/A
Energy consumption - default setpoint	1.5 kWh/24h
Heat emission 100%	198 Watt
Heat emission default setpoint	N/A
Nominal consumption	288 Watt/1.4 A
Start current	10.1 A
Sound level	45.5 dB(A)



BIOBASIC RF600

Technical specifications	Data
Temperature range	-25°C/-5°C
SW variant	F71
Defrost system	Automatic smart defrost with re-evaporation of the defrost water. Heating element
ATEX certificate	DTI 19ATEX0114X
ATEX Marking	II 3G Ex ec nC ic IIB T1 Gx
Refrigerant	R290
Refrigerant charge	73 g
Refrigeration capacity at -10°C	339 Watt
GWP – CO2e	N/A
Energy consumption – default setpoint	3.9 kWh/24h
Heat emission 100%	344 Watt
Heat emission default setpoint	N/A
Nominal consumption	357 Watt/1.8 A
Start current	13.8 A
Sound level	44.3 dB(A)

Declaration of Conformity



English EC Declaration of Conformity

We, **Gram Commercial** declare under sole responsibility that the following products:

Range:	BIOBASIC
Model:	210, 310, 410 & 600
Refrigeration:	R600a, R290 & R134a
Product description:	Refrigerators and freezer for laboratory
Valid from (Year/Week):	2022/02

To which this declaration relates, is in compliance with all the applicable essential requirements, and other provisions of the European Council Directive and regulations.

Directives and Regulations of the European Parliament and of the Council:

Machinery Directive 2006/42/EC
 - ATEX Directive 2014/34/EU
 - Pressure Equipment Directive 2014/68/EU
 - Low Voltage Directive 2014/35/EU
 - EMC Directive 2014/30/EU
 - RoHS Directive 2011/65/EU
 - REACH EC No.1907/2006
 - F-Gas Regulation (EU) No 517/2014

Product compliance has been demonstrated on the basis of:

Harmonized Standards:	Text:
EN 61010-1:2010	Safety requirements for electrical equipment for measurement, control, and laboratory use - Part 1: General requirements
EN 61326-1:2013	Electrical equipment for measurement, control and laboratory use – EMC requirements – Part 1: General requirements
EN IEC 60079-0:2018 EN IEC 60079-0:2018/AC:2020	Explosive atmospheres – Part 0: Equipment – General requirements
EN 60079-7:2015 EN 60079-7:2015/A1:2018	Explosive atmospheres – Part 7: Equipment protection by increased safety "e"
EN 60079-11:2012	Explosive atmospheres – Part 11: Equipment protection by intrinsic safety "i"
EN IEC 60079-15:2019	Explosive atmospheres – Part 15: Equipment protection by type of protection "n"
EN 60079-18:2015	Explosive atmospheres – Part 18: Equipment protection by encapsulation "m"
EN ISO 3744:2010	Acoustics – Determination of sound power levels and sound energy levels of noise sources using sound pressure – Engineering methods for an essentially free field over a reflecting plane
EN ISO 9001:2015	Quality management systems – Requirements
EN ISO 14001:2015	Environmental management systems – Requirements with guidance for use

Gram Commercial
 Aage Grams Vej 1
 DK-6500 Vojsens
 Telephone: + 45 73 20 13 00


Vojsens, 11.1.2022

John B. S. Petersen
 Approval Manager


Rev. 005 – 11.1.2022

Wiring diagrams


BIOBASIC RR/RF210, RR/RF310, RR/RF410

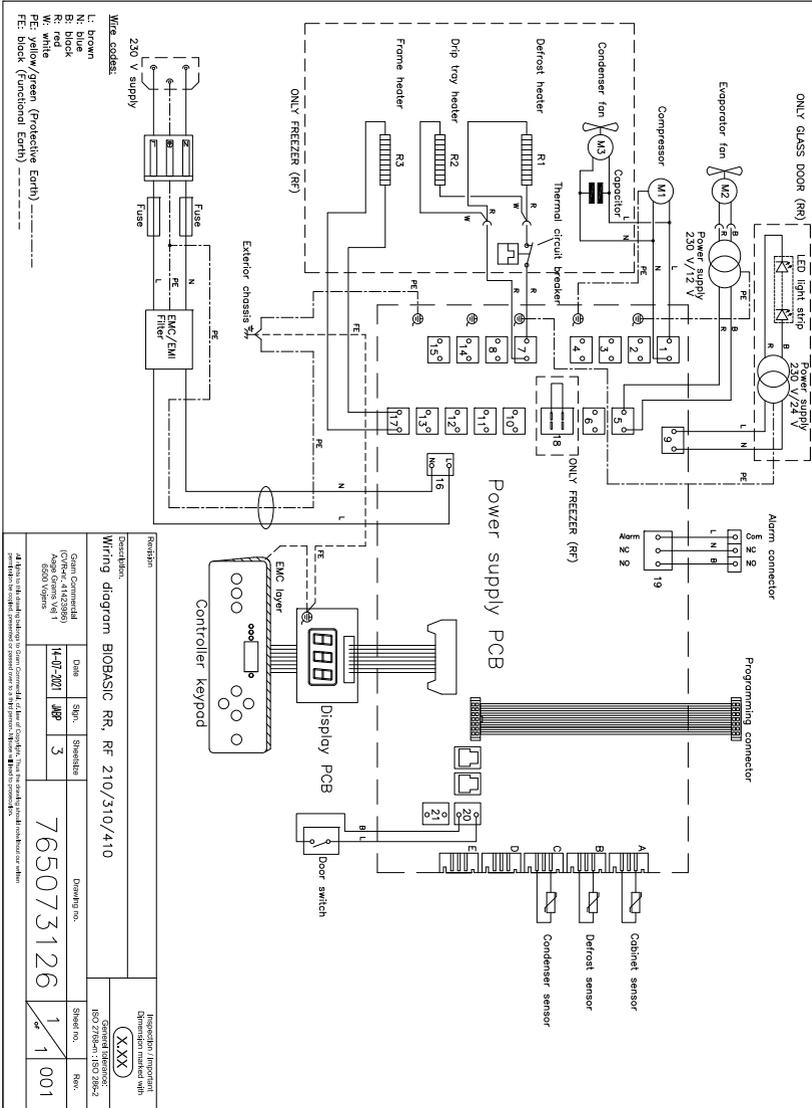


BIOBASIC RR600

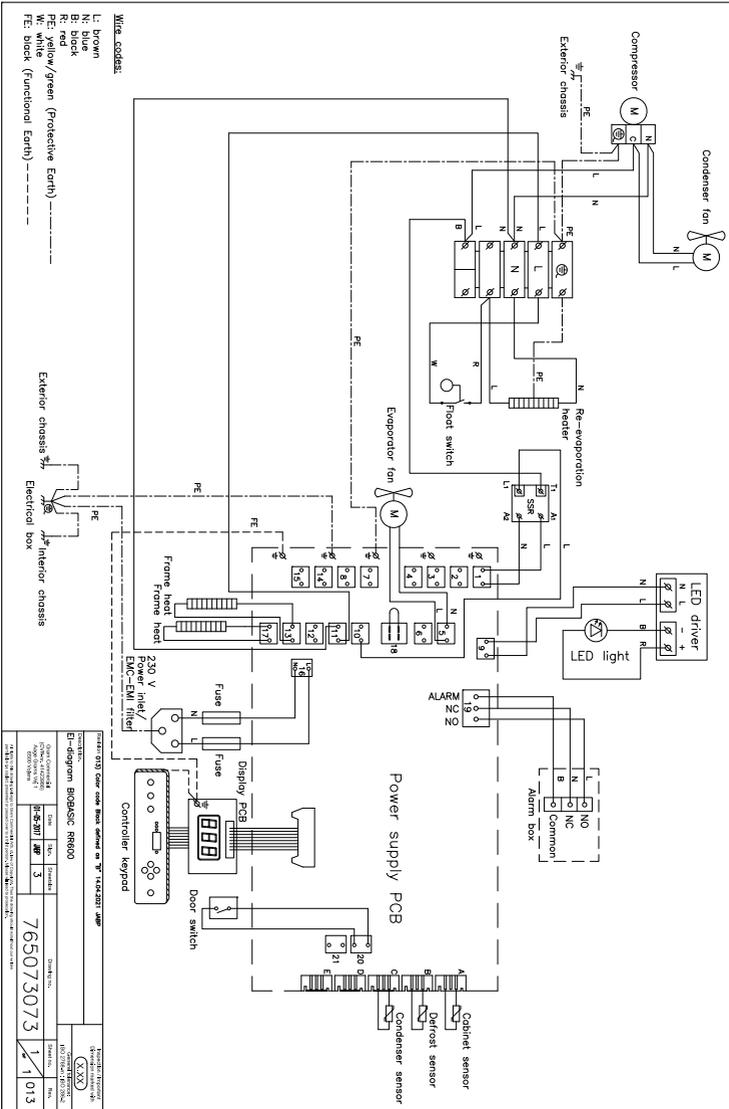

BIOBASIC RF600

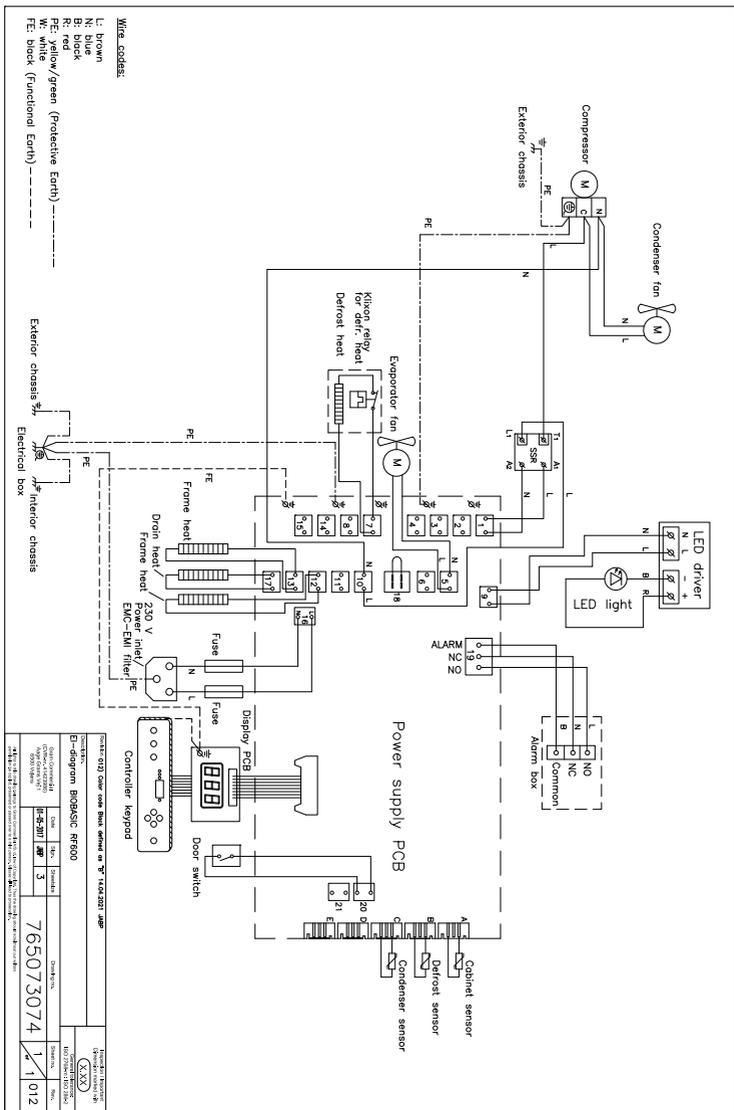

Piping diagram

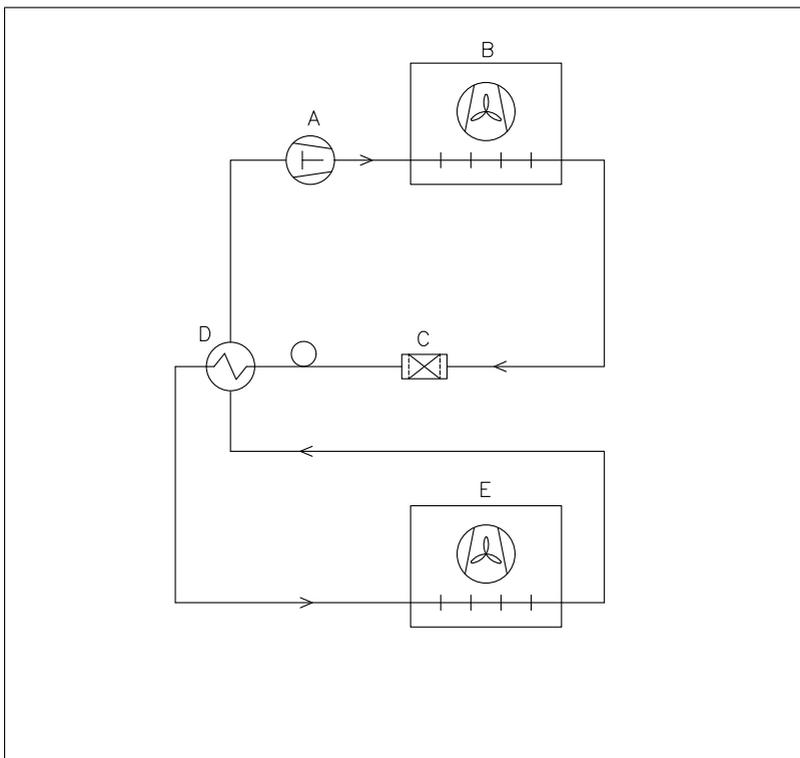

	DK	GB	D
A	Kompressor	Compressor	Kompressor
B	Kondensator	Condenser	Verflüssiger
C	Tørrefilter	Filter drier	Trockenfilter
D	Varmeudveksler	Heat exchanger	Wärmeaustauscher
E	Fordamper	Evaporator	Verdampfer
Revision			Inspection / important Dimension marked with
Description: Piping diagram BioCompact II			(X.XX)
			General tolerance: ISO 2768-m : ISO 286-2
Gram Commercial (CVR-nr. 41423986) Aage Grams Vej 1 6500 Vojens	Date	Sign.	Sheetsize
	18-01-2022	JABP	4
Drawing no. 765042593			
		Sheet no. 1	Rev. 000
		of 1	
All rights to this drawing belongs to Gram Commercial A/S. of. Inve of Copyright. Thus the drawing should notwithout our written permission be copied, presented or passed over to a third person. Misuse will lead to prosecution.			

Notes





Manufactured by Gram Commercial

Aage Grams Vej 1 · 6500 Vojens · Denmark

Tel: +45 73 20 13 00

e-mail: info@gram-bioline.com · www.gram-biobasic.com

Gram Commercial, filial af Hoshizaki Europe BV, Holland



Item number **765042590** Rev.: **20220926_1**

Language: **English**