

Spare parts – BioCompact II
RR 610 H
Glass door




Description :

**Spare parts list for BioCompact II
RR 610 H Glass door**

Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
25.01.12	JP	19.09.25	jolu	016		

Exploded view

0925 -

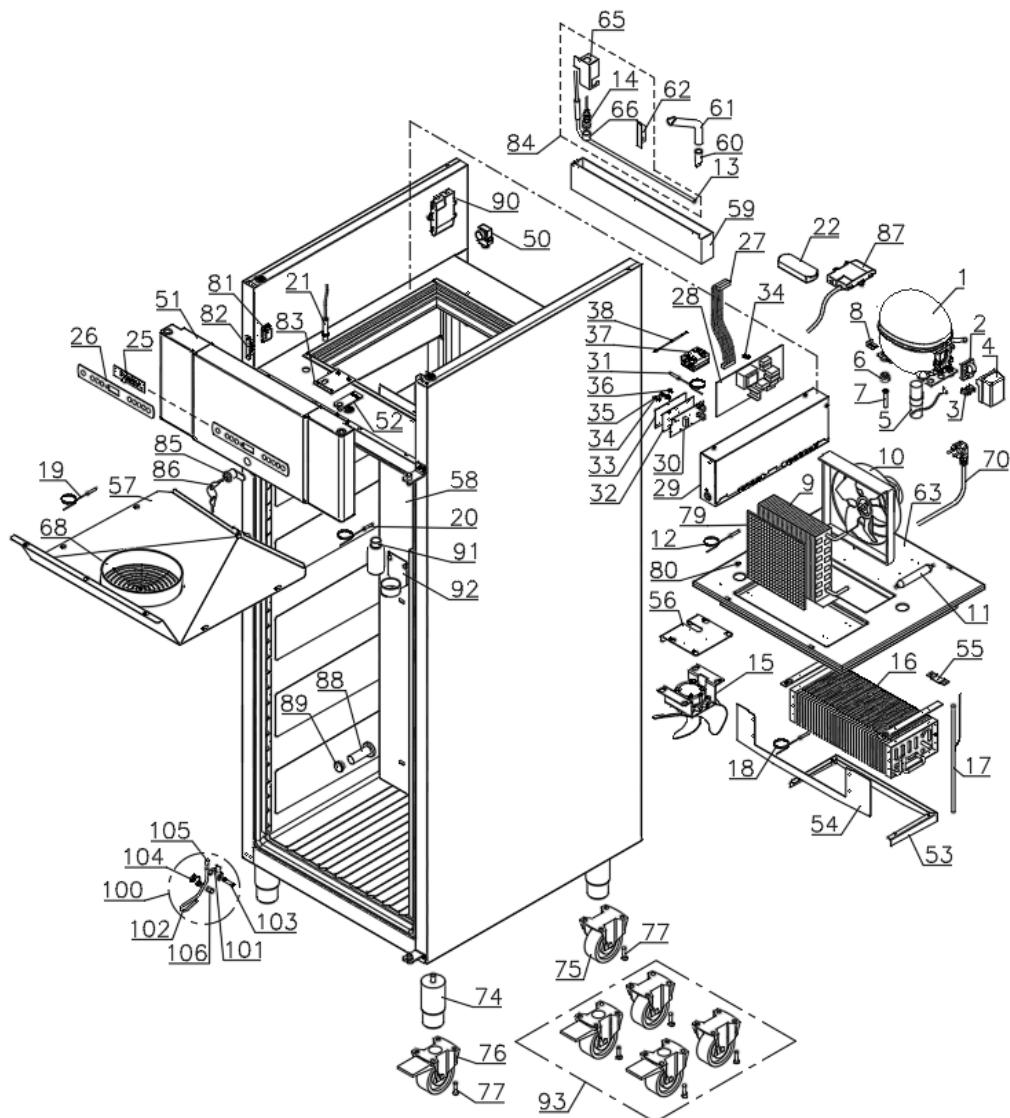


Description :

Spare parts list for BioCompact II
RR 610 H Glass door

Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
25.01.12	JP	19.09.25	jolu	016		

Glass door

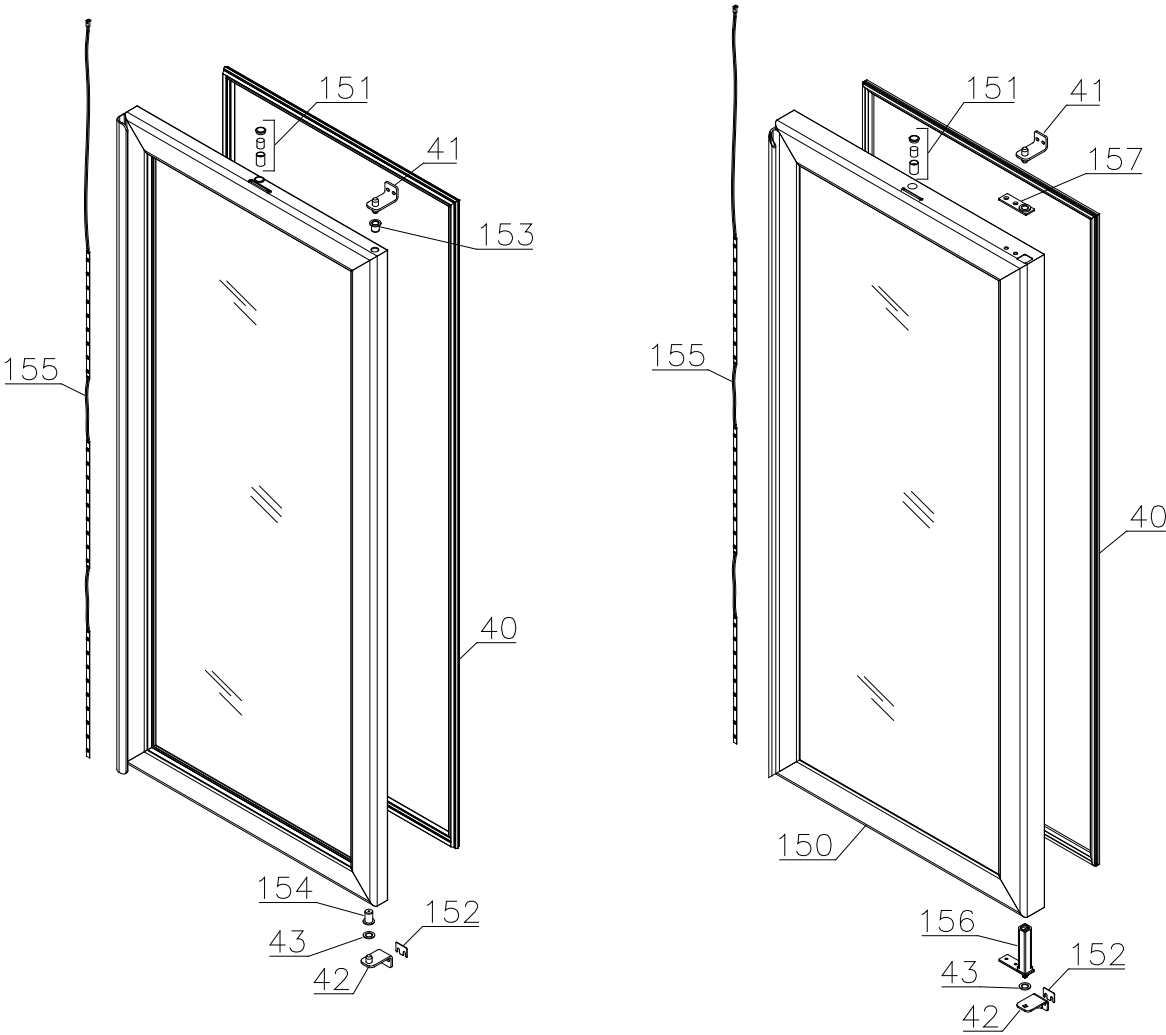


Description :

**Spare parts list for BioCompact II
RR 610 H Glass door**

Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
25.01.12	JP	19.09.25	jolu	016		

Interior parts

Shelves:

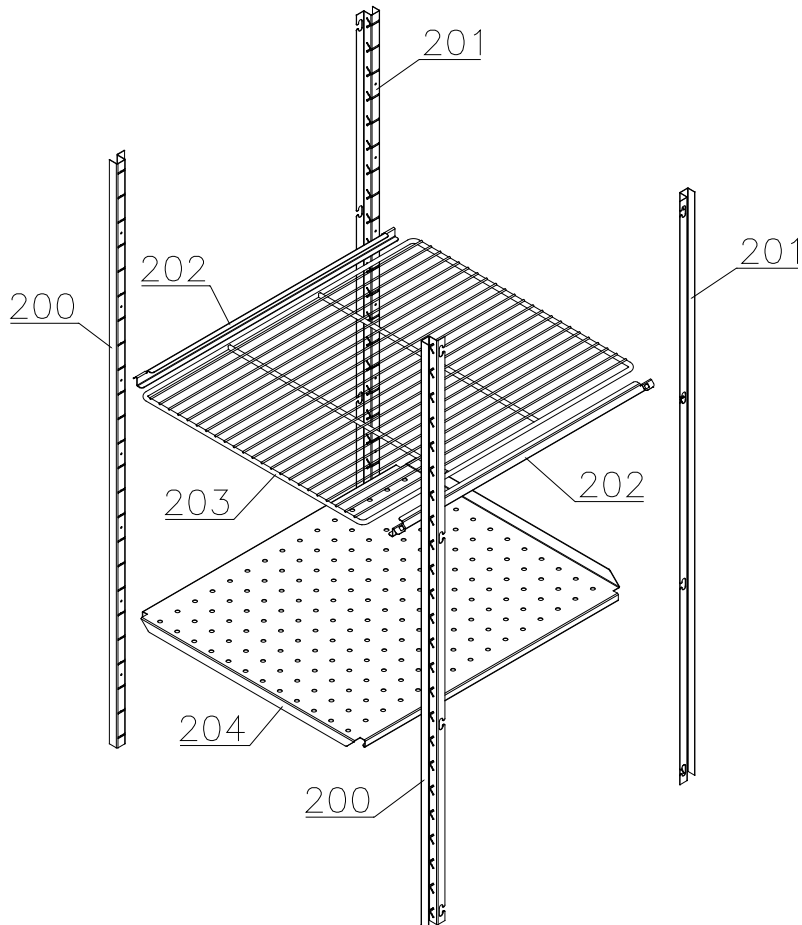


Description :

**Spare parts list for BioCompact II
RR 610 H Glass door**

Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
25.01.12	JP	19.09.25	jolu	016		

Drawers:

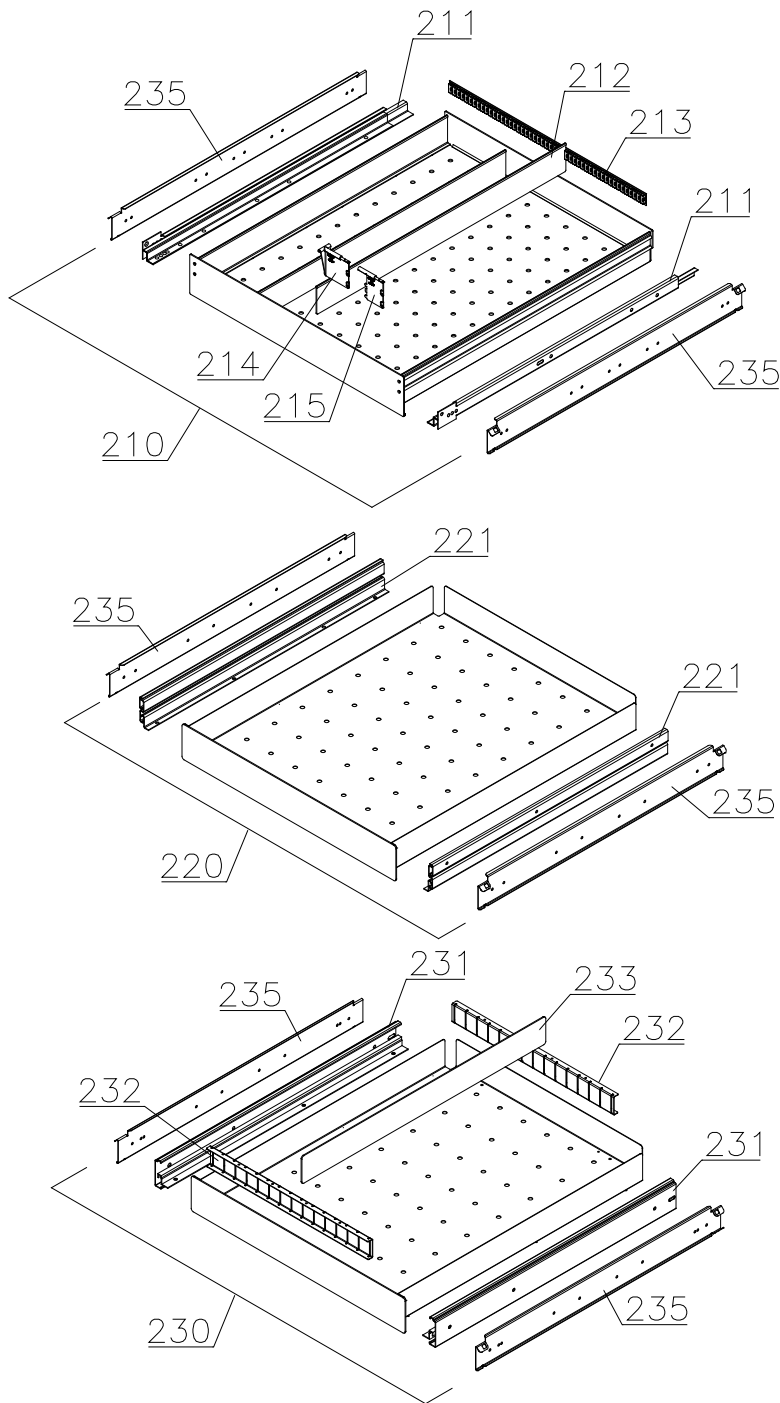

Gram Commercial Technical information					765041486
Description : Spare parts list for BioCompact II RR 610 H Glass door					
Compiled.		Latest revision.			Deleted.
Date	Name	Date	Name	Rev.no	Date
25.01.12	JP	19.09.25	jolu	016	

Item no.	Description	Gram part no.	week/year introduced	week/year deleted
----------	-------------	---------------	----------------------	-------------------

Essential spare parts

1	Compressor FR 7.5 G – 50/60 Hz	500550061	1508	
2	Starting relay	410589015	1508	
3	Cord relief	452440297	1508	
4	Cover	452440293	1508	
5	Start capacitor	451720018	1508	
6	Rubber foot	760080519	1508	
7	Screw for compressor	423200200	1508	
8	Fastening clip	760082367	1508	
9	Condenser	612900500	1508	
10	Condenser fan guard compl.	765080189	1508	
11	Filter drier	762900157	1508	
12	Condenser sensor	453531500	1508	
13	Re-evaporation heater	765071848	1508	2714
13	Re-evaporation heater	765072862	2714	
14	Float switch	765072049	1508	
15	Evaporator fan complete	761221902	1508	
16	Evaporator	761220332	1508	
17	Heat exchanger	761221403	1508	
18	Defrost sensor	453531501	1508	
19	Cabinet sensor	453531501	1508	
20	E-sensor	453531504	1508	
21	Door switch	450400113	1508	
22	LED driver	765072844	1508	
25	Controller display part	612900056	0508	0825
25	Controller display part	612900059	0525	
26	Label with keys	765041753	0508	0825
26	Label with keys	765041889	0925	
27	Band cable	453582046	1508	
27	Band cable for LTP	765072958	3519	
28	Electronic controller K6+	761256154	0508	
29	Electrical box complete	761232111	0508	0825
29	Electrical box complete	761233701	0925	
29	Electrical box complete (LTP)	761233703	0925	
30	LTP controller	453580515	3519	
30	EMC/LTP kit - retro fit BioCompact 610	761234073	3519	0925
30	EMC/LTP kit - retro fit BioCompact 610	761234089	1025	
31	LTP sensor	453531526	3519	
32	PCB protection sheet	765072698	0718	
33	Screening plate	765073271	0718	
34	Dual locking spacer H=7.9 mm	451119525	0718	0825
34	Dual locking spacer H=5.0 mm	451119520	0925	
35	Dual locking spacer H=15.9 mm	451119526	0718	
36	Push spacer H=7.9 mm	451119527	0718	0825
36	Push spacer H=5.0 mm	451119533	0925	

week/year week/year

Gram Commercial Technical information					765041486
Description : Spare parts list for BioCompact II RR 610 H Glass door					
Compiled.		Latest revision.			Deleted.
Date	Name	Date	Name	Rev.no	Date
25.01.12	JP	19.09.25	jolu	016	

Item no.	Description	Gram part no.	introduced	deleted
37	Solid state relay	450200300	4717	
38	Cable to SSR incl. resistance	765071445	1023	
38	Cable to SSR incl. resistance "LTP"	765071449	1023	
40	Door gasket	761220372	1508	4314
40	Door gasket	761222871	4314	
41	Hinge, top	761222735	1508	
42	Hinge, bottom	761222737	1508	4314
42	Hinge, bottom self-close right	761222824	4314	
42	Hinge, bottom self-close left	761222827	4314	
43	Washer	760250524	1508	4314
43	Washer	420178003	4314	

Gram Commercial Technical information					765041486
Description : Spare parts list for BioCompact II RR 610 H Glass door					
Compiled.		Latest revision.			Deleted.
Date	Name	Date	Name	Rev.no	Date
25.01.12	JP	19.09.25	jolu	016	

Item no.	Description	Gram part no.	week/year introduced	week/year deleted
----------	-------------	---------------	----------------------	-------------------

Other spare parts

50	Air relief valve	765081206	1508	
51	Front panel	stainless 761234020	0508	0825
51	Front panel	stainless 761234020	0925	
51	Front panel	white 761234021	0508	0825
51	Front panel	white 761234021	0925	
51	Front panel Chart Recorder	white 761234169	0925	
51	Front panel Chart Recorder	stainless 761234170	0925	
52	Hinge for front panel	761220388	1508	
53	Evaporator cover, side/back	761220326	1508	
54	Evaporator cover, front	761220325	1508	
55	Suspension bracket	760250295	1508	
56	Fan holder	761220369	1508	
57	Drip water tray	761221430	1508	
58	Air distribution plate	761220331	1508	
59	Re-evaporation tray	765081149	1508	
60	Water trap	760082891	1508	
61	Drain pipe	765081153	1508	5017
61	Drain pipe	765081253	5017	
62	Heater bracket	761255445	1508	2714
62	Heater bracket	763250002	2714	
63	Insulation slab	761221734	1508	
64	Screw for fan blade	612900100	1508	
65	Float switch bracket	761255446	1508	2714
65	Float switch bracket	763250003	2714	
66	Plastic cap	450450080	1508	
68	Fan guard	765081239	1508	
70	Lead and plug, DK 3-pin	765072568	1508	
70	Lead and plug, Schuko	765072566	1508	
70	Lead and plug, Switzerland	765070163	1508	
70	Lead and plug, GB	765072567	1508	
74	Leg, stainless H=135-200	765080115	1508	
75	Fixed wheel, H=125	480270520	1508	
76	Castor wheel, with brake	480260050	1508	
77	Screw M8 x 12	422715000	1508	
79	Condenser filter	761114416	1508	
80	Attachment plug	765081246	2809	
81	Front panel close catch	761220375	1508	
82	Front panel close keeper	761220374	1508	
83	Bracket for lock	761220376	1508	
84	Re-evap. heater kit complete	761255438	1508	2714
84	Re-evap. heater kit complete	763254026	2714	

Gram Commercial Technical information					765041486
Description : Spare parts list for BioCompact II RR 610 H Glass door					
Compiled.		Latest revision.			Deleted.
Date	Name	Date	Name	Rev.no	Date
25.01.12	JP	19.09.25	jolu	016	

Item no.	Description	Gram part no.	week/year introduced	week/year deleted
85	Door lock	765080092	1508	
86	Key no. 1 (25010)	765080340	1508	
86	Key no. 2 (25020)	765080341	1508	
86	Key no. 3 (25030)	765080342	1508	
86	Key no. 4 (25040)	765080343	1508	
86	Key no. 5 (25050)	765080344	1508	
86	Master key	765080385	1508	
87	Box for alarm connection	761233402	0508	
88	Insulation plug	765080563	0508	
89	Plastic cap	501020023	0508	
90	Electrics cover	765070143	1508	
91	Liquid container	765080384	1508	
92	Holder for liquid container	761256497	1508	
93	Set of castor wheels	818899905	1508	
100	Foot pedal complete	819581053	0911	
101	Bracket for pedal	761224010	0911	
102	Pedal arm	761224002	0911	
103	Spindle	761224007	0911	
104	Spring	761224005	0911	
105	Plastic cap	581400021	0911	
106	Bush	761224008	0911	
150	Glass door, self-close	761232965	1508	
151	Door magnet complete	761250786	1508	
152	Spacer f. hinge	761258621	1508	
153	Bush, top	760250525	1508	4314
154	Bush, bottom	761220936	1508	4314
155	LED light strip	761222718	1508	
156	Self-close device	763251745	4314	
157	Hinge bearing complete	761220980	4314	

Gram Commercial Technical information					765041486
Description : Spare parts list for BioCompact II RR 610 H Glass door					
Compiled.		Latest revision.			Deleted.
Date	Name	Date	Name	Rev.no	Date
25.01.12	JP	19.09.25	jolu	016	

Item no.	Description	Gram part no.	week/year introduced	week/year deleted
----------	-------------	---------------	----------------------	-------------------

Interior equipment:

200	Wall rail, front	761220335	1508	
201	Wall rail, back	761220336	1508	
202	Shelf support rail	761256455	1508	
203	Grey wire shelf	818880008	1508	
204	Perforated shelf, stainless	761230083	0508	
210	Alu drawer, complete	761230362	2210	
211	Set of sliding rails	761234055	2210	
212	Divider rail	761230016	2210	
213	Holder for divider rail	761230063	2210	
214	Plastic divider 1	761230021	2210	
215	Plastic divider 2	761230022	2210	
220	Stainless drawer 30 kg, complete	761231232	1508	
221	Set of sliding rails	761234000	1508	
230	Stainless drawer 40 kg, complete	761231202	1508	
231	Set of sliding rails	761234058	1508	
232	Divider rail	761230305	1508	
233	Divider	761230308	1508	
235	Support for sliding rail	761230201	1508	3310
235	Support for sliding rail, right	761230278	3310	
235	Support for sliding rail, left	761230281	3310	