

Spare parts

BioUltra

UL 570




Description :

**Spare parts list
BioUltra UL 570**

Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
05.07.13	JP	01.12.17	JP	006		

Compressor unit

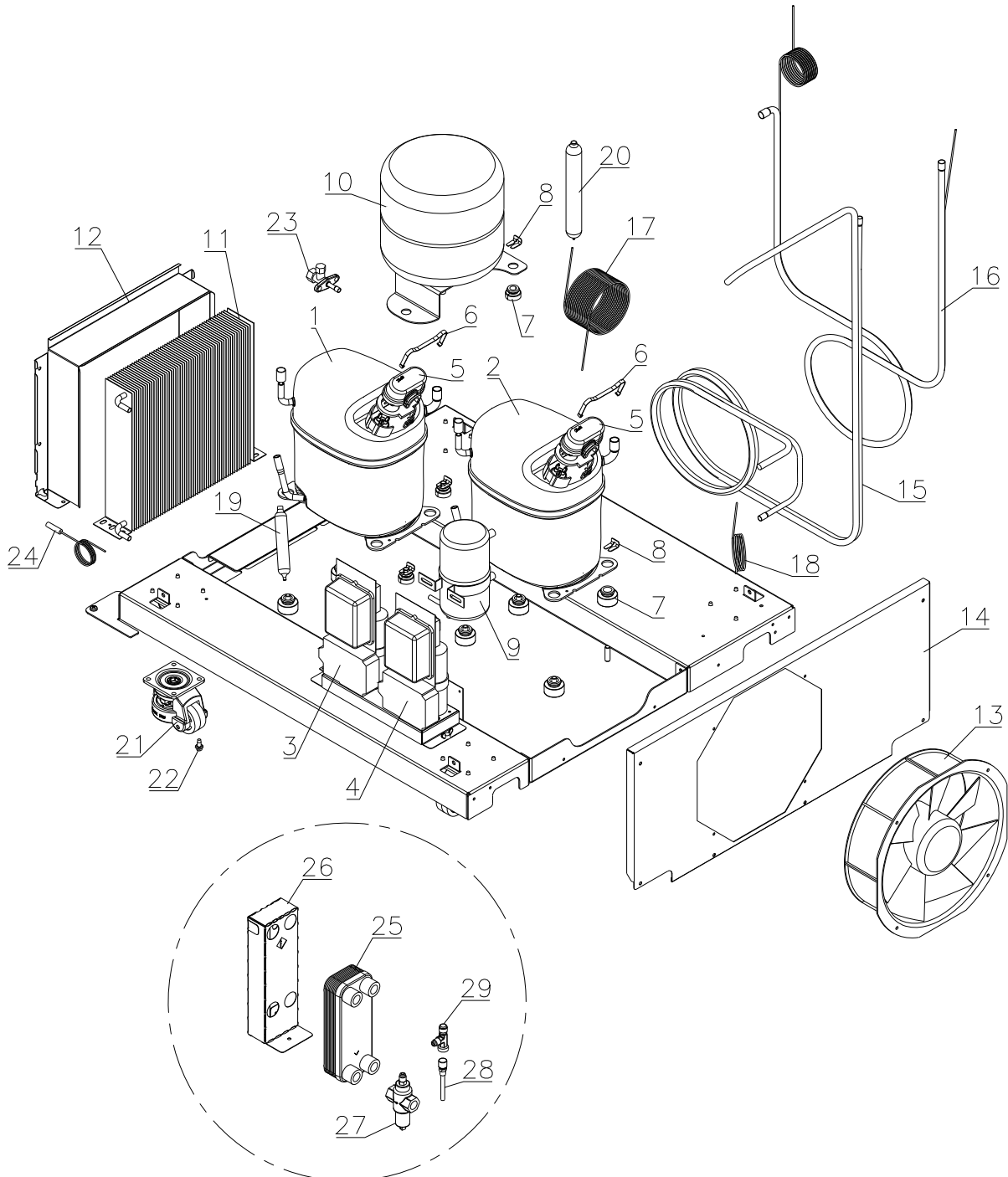


Description :

**Spare parts list
BioUltra UL 570**

Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
05.07.13	JP	01.12.17	JP	006		

Electrical parts

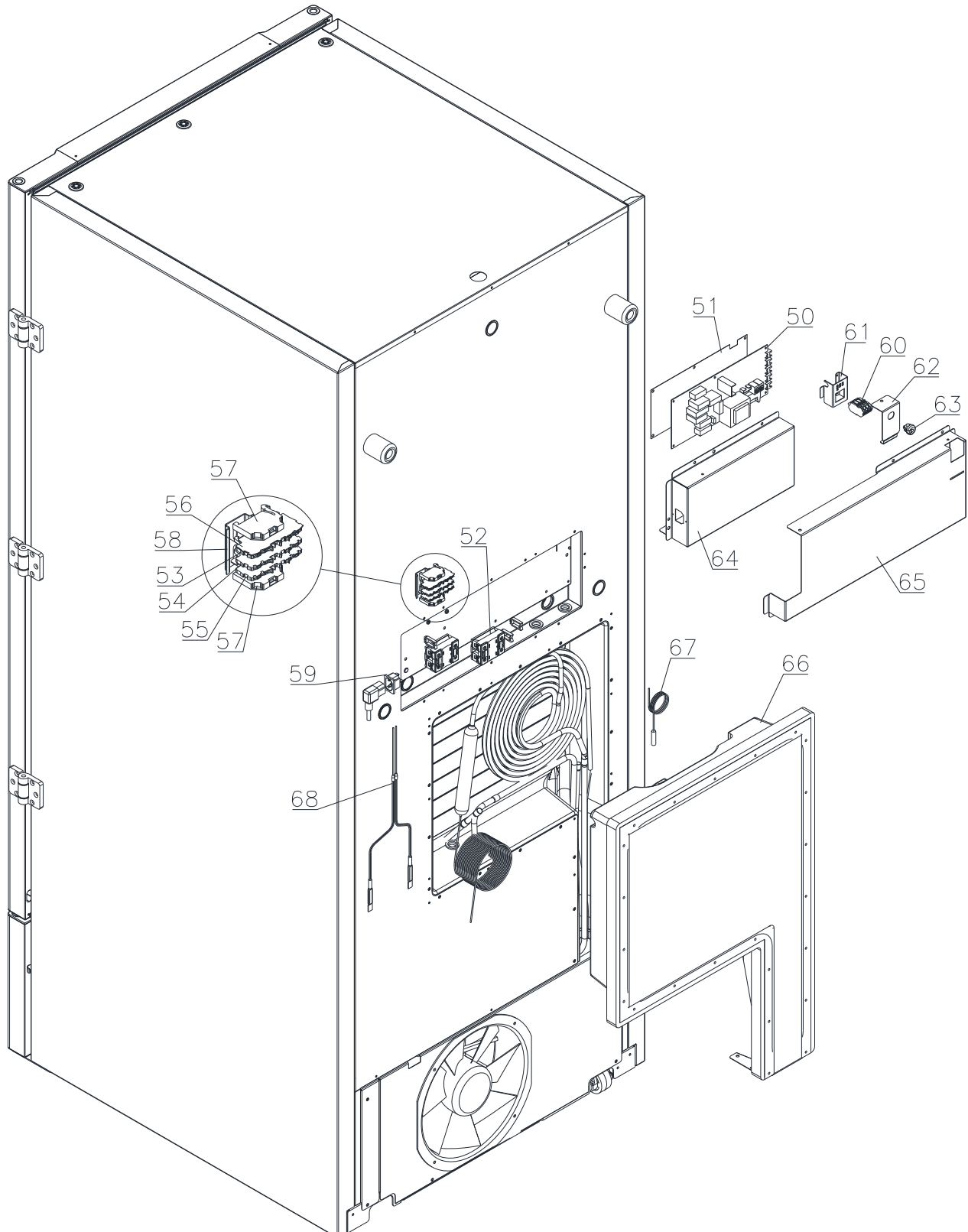


Description :

**Spare parts list
BioUltra UL 570**

Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
05.07.13	JP	01.12.17	JP	006		

Cabinet parts

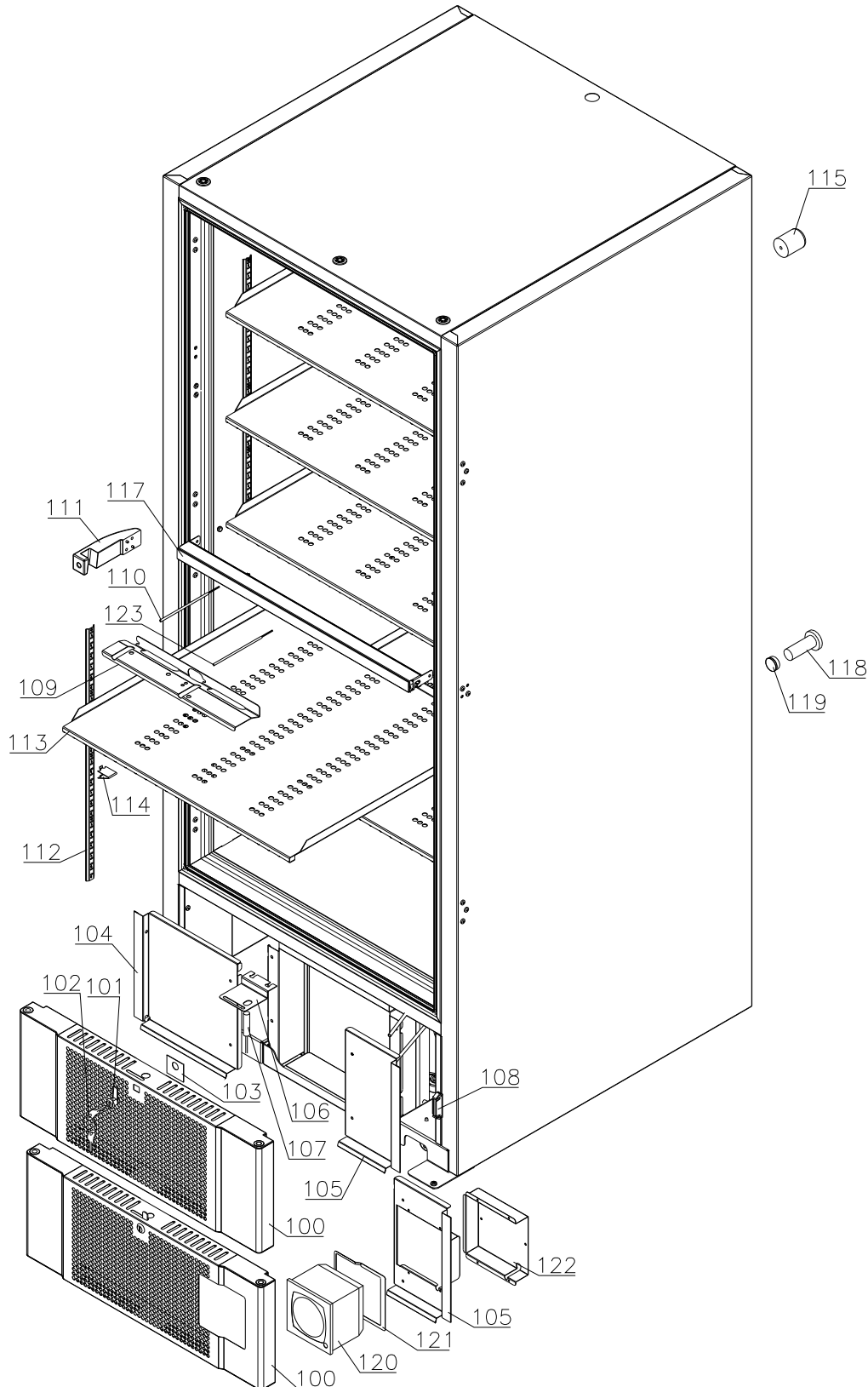


Description :

**Spare parts list
BioUltra UL 570**

Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
05.07.13	JP	01.12.17	JP	006		

Internal doors

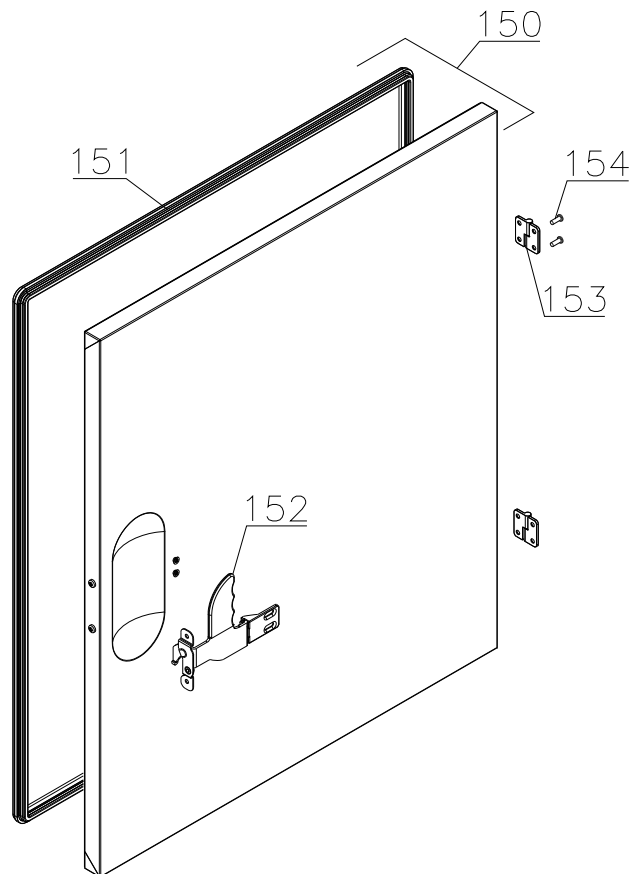


Description :

Spare parts list
BioUltra UL 570

Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
05.07.13	JP	01.12.17	JP	006		

External door

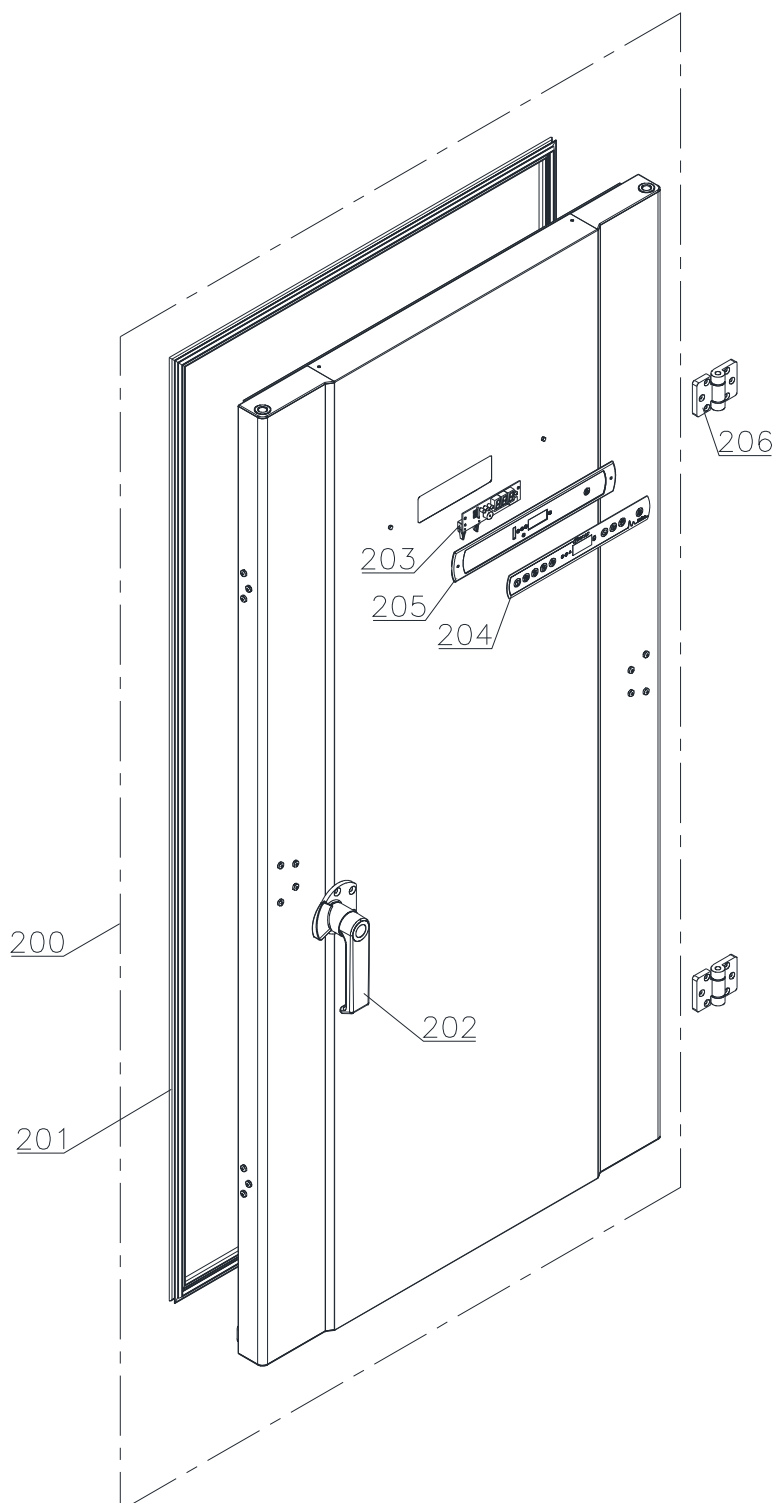

Gram Commercial Technical information					765041607	
<u>Description :</u> Spare parts list BioUltra UL 570						
Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
05.07.13	JP	01.12.17	JP	006		

Item no.	Description	Gram part no.	W/Y Issued	Deleted
<u>Compressor unit</u>				
1	Compressor GS26CLX (HT)	500550600	0114	
2	Compressor GS34CLX (LT)	500550601	0114	2115
2	Compressor GS26CLX (LT)	500550600	2115	
3	Starting device (GS26CLX)	453051005	0114	
4	Starting device (GS34CLX)	453051006	0114	2115
5	Cover	452440281	0114	
6	Clamp	452440284	0114	
7	Rubber bush	760080519	0114	
8	Compressor clip	760082367	0114	
9	Oil separator	600020052	0114	
10	Expansion tank	600020020	0114	
11	Condenser	765080166	0114	
12	Condenser shroud	761238666	0114	
13	Condenser fan	761238589	0114	
14	Fan shroud	761238669	0114	
15	Heat exchanger LT	760000337	0114	
16	Heat exchanger HT	760000338	0114	
17	Capillary tube LT	760000341	0114	
18	Capillary tube L=700 mm	760000340	0114	0916
19	Filter drier HT	765080158	0114	
20	Filter drier LT	765080194	0114	
21	Castor, adjustable	480260105	0114	
22	Screw M6x12	422712010	0114	
23	Auxiliary valve	460978041	1716	
24	Condenser sensor	453531504	0114	
<u>Water cooled condenser parts</u>				
25	Condenser 2 kW	765080221	1817	
26	Condenser holder	761238588	1817	
27	Water valve	765080149	1817	
28	Schrader valve	765080148	1817	
29	T-nipple w/schrader valve	765080519	1817	

Gram Commercial Technical information					765041607	
<u>Description :</u> Spare parts list BioUltra UL 570						
Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
05.07.13	JP	01.12.17	JP	006		

Item no.	Description	Gram part no.	W/Y Issued	Deleted
<u>Electrical parts / cascade system</u>				
50	Controller power part	761256142	0114	
51	Plastic foil	765040777	0114	
52	Contactor B6-40-00	450241000	0114	3515
52	SSR relay	450200300	3515	
53	Terminal block, grey	765070706	0114	
54	Terminal block, blue	765070707	0114	
55	Terminal block, yellow/green	765070708	0114	
56	Terminal block end plate	765070709	0114	
57	Terminal block end stop	765070711	0114	
58	DIN rail	760254475	0114	
59	Appliance inlet	451022002	0114	
60	Terminal block for alarm	765070705	0114	
61	Bracket for alarm terminal	761232157	0114	
62	Cover for alarm terminal	761238673	0114	
63	Cord relief	422210002	0114	
64	Cover for high current parts	761238672	0114	
65	Cover for low current parts	761238677	0114	
66	Cover for cascade system	761238682	0114	
67	Cascade sensor	453531505	0114	
68	Defrost heater	765072758	0114	

Gram Commercial Technical information					765041607	
<u>Description :</u> Spare parts list BioUltra UL 570						
Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
05.07.13	JP	01.12.17	JP	006		

Item no.	Description	Gram part no.	W/Y Issued	Deleted
<u>Cabinet parts</u>				
100	Front panel complete, white	761231490	0114	
100	Front panel complete, stainless	761234033	0114	
100	Front panel compl, f. chart rec., white	761234034	0114	
101	Lock	765080092	0114	
102	Key no. 1 (25010)	765080340	0114	
102	Key no. 2 (25020)	765080341	0114	
102	Key no. 3 (25030)	765080342	0114	
102	Key no. 4 (25040)	765080343	0114	
102	Key no. 5 (25050)	765080344	0114	
102	Master key	765080385	0114	
103	Lock bracket	760249008	0114	
104	Condenser screen, left	761238664	0114	
105	Condenser screen, right	761238663	0114	
105	Condenser screen, right, f. ch.recorder	761238571	0114	
106	Bracket for door switch	761238562	0114	
107	Door switch	450400102	0114	
108	Magnet for front panel	761253312	0114	
109	Sensor bracket	761238553	0114	
110	Room sensor	453531513	0114	
111	Strike side-mount, righthand	765080471	0114	
111	Strike side-mount, lefthand	765080473	0114	
112	Wall rail	761238594	0114	
113	Shelf	761238599	0114	
114	Shelf support	761238592	0114	
115	Spacer bush L=50	762251184	0114	
117	Cabinet crossbar	761238628	0114	
118	Insulation plug	765080563	0114	
119	Plastic cap	501020023	0114	
120	Chart recorder	761256515	0114	
121	Chart recorder bracket	761238569	0114	
122	Back cover	761238570	0114	
123	E-sensor	453531513	0114	

Gram Commercial Technical information					765041607	
<u>Description :</u> Spare parts list BioUltra UL 570						
Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
05.07.13	JP	01.12.17	JP	006		

Item no.	Description	Gram part no.	W/Y Issued	Deleted
<u>Internal door parts</u>				
150	Internal door complete	761234019	0114	
151	Gasket	761238649	0114	
152	Handle	761238555	0114	
153	Hinge	424180032	0114	
154	Screw	420182025	0114	
<u>External door parts</u>				
200	Door complete, white	761231489	0114	
200	Door complete, stainless	761234035	0114	
201	Gasket	761238641	0114	
202	Handle, righthand	765080470	0114	
202	Handle, lefthand	765080472	0114	
203	Electronic control, display part	612900052	0114	
204	Label with keys	765041752	0114	
205	Display holder	761238685	0114	
206	Hinge	424180030	0114	