

Spare parts – BioCompact II

RR 610 G




Spare parts list for BioCompact II RR 610 G

Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
09.01.09	JP	19.09.25	Jolu	027		

Exploded view RR 610

0508 – 0825

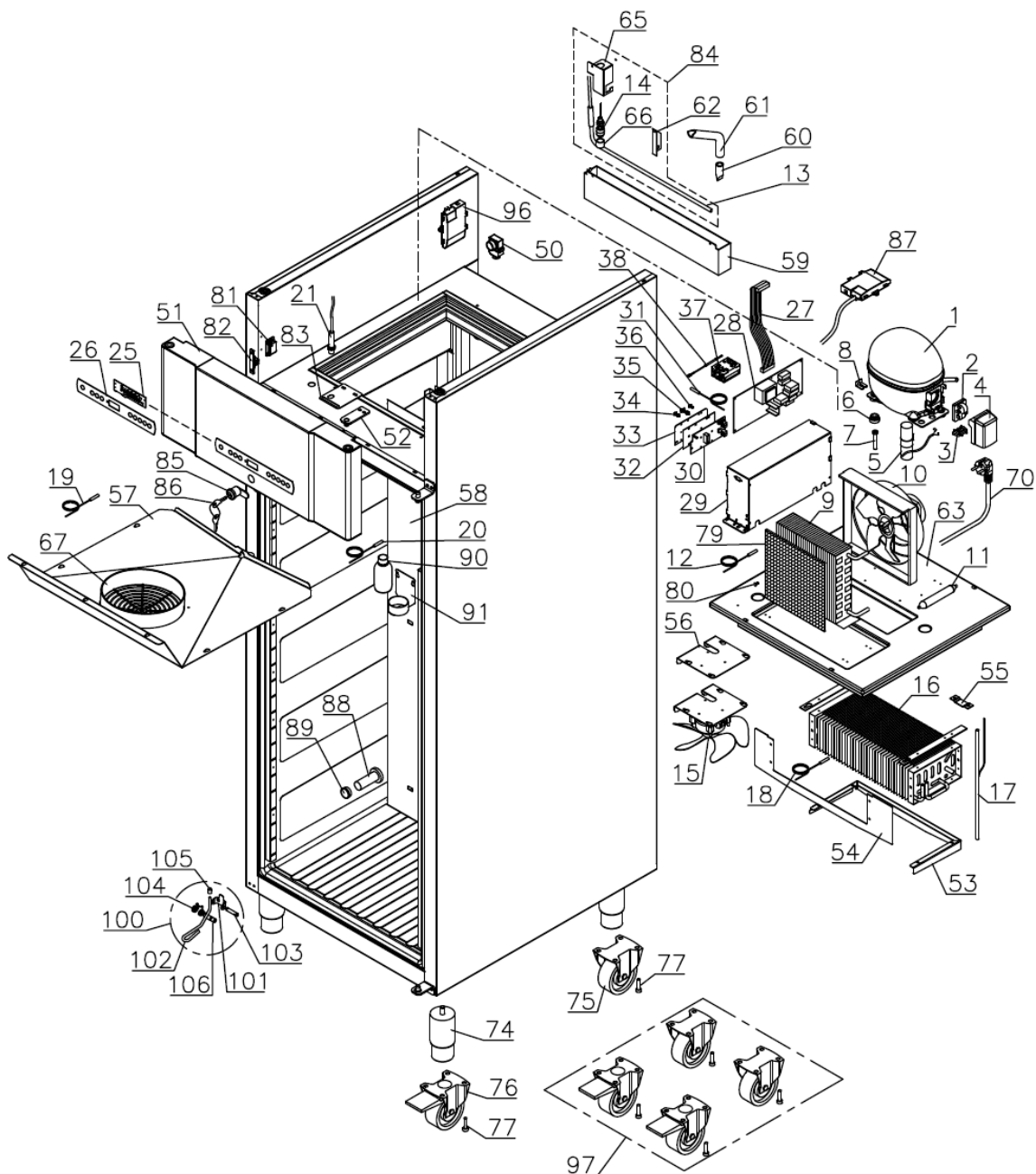


Description :

**Spare parts list for BioCompact II
RR 610 G**

Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
09.01.09	JP	19.09.25	Jolu	027		

Exploded view RR 610


0925 -

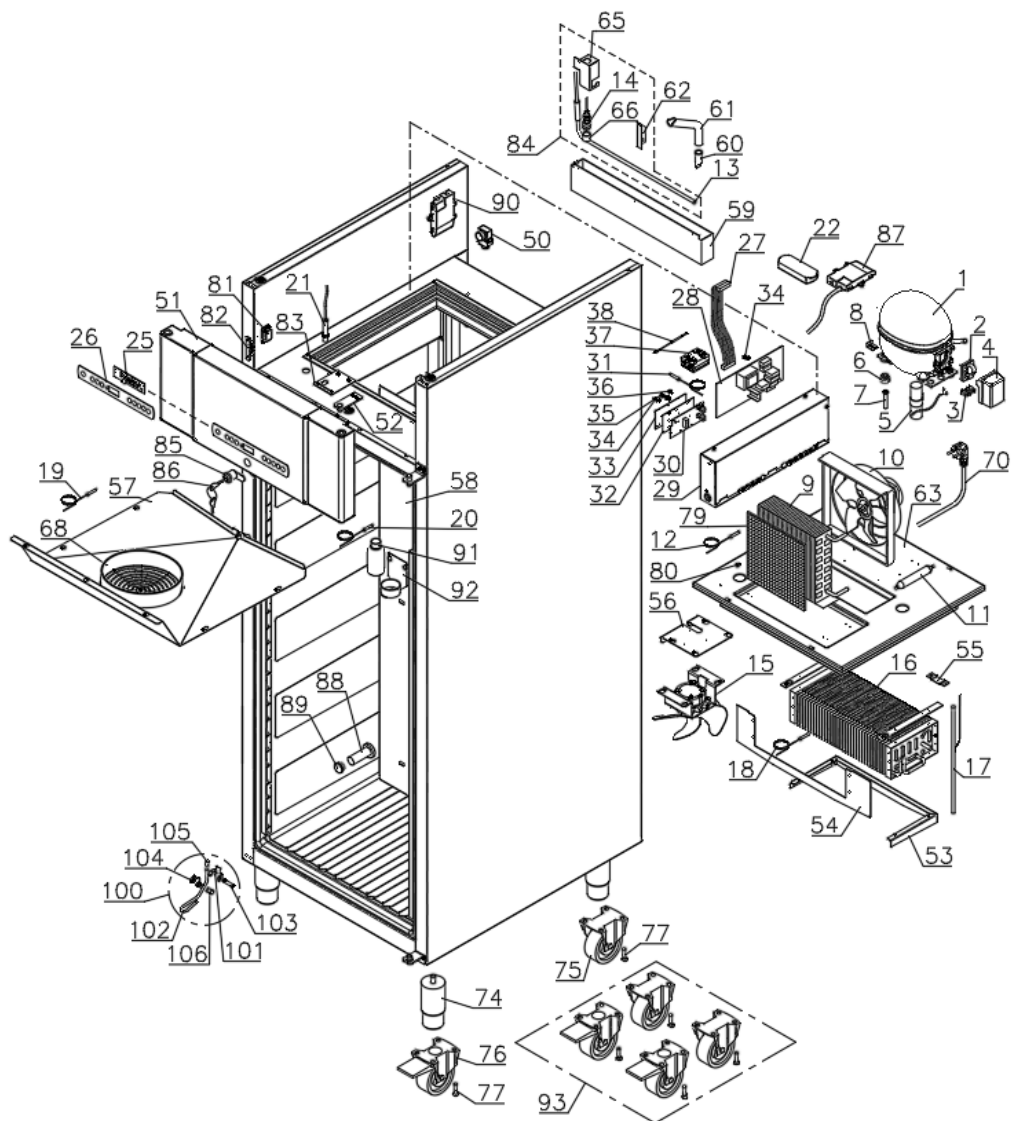

Technical information					765041294	
<u>Description :</u> Spare parts list for BioCompact II RR 610 G						
Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
09.01.09	JP	19.09.25	Jolu	027		


Door

0508-4314

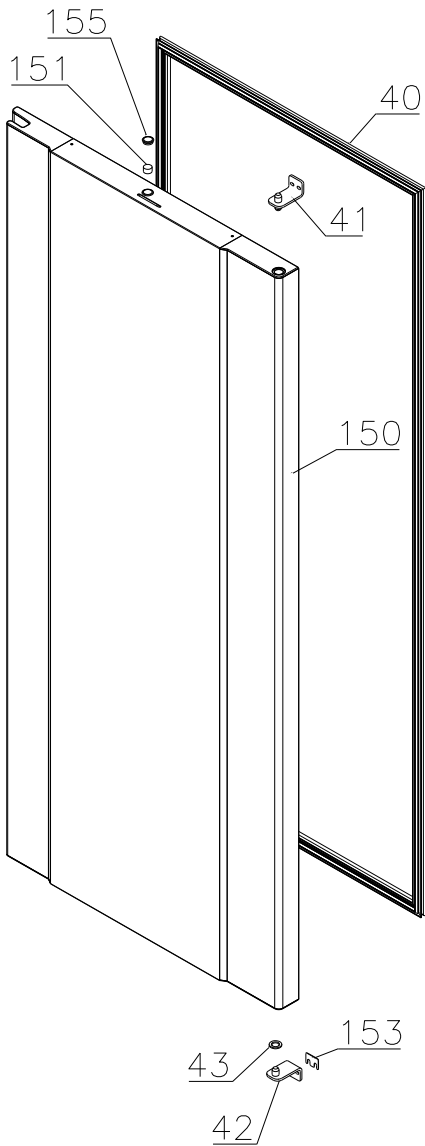

Technical information					765041294	
<u>Description :</u> Spare parts list for BioCompact II RR 610 G						
Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
09.01.09	JP	19.09.25	Jolu	027		

4314-

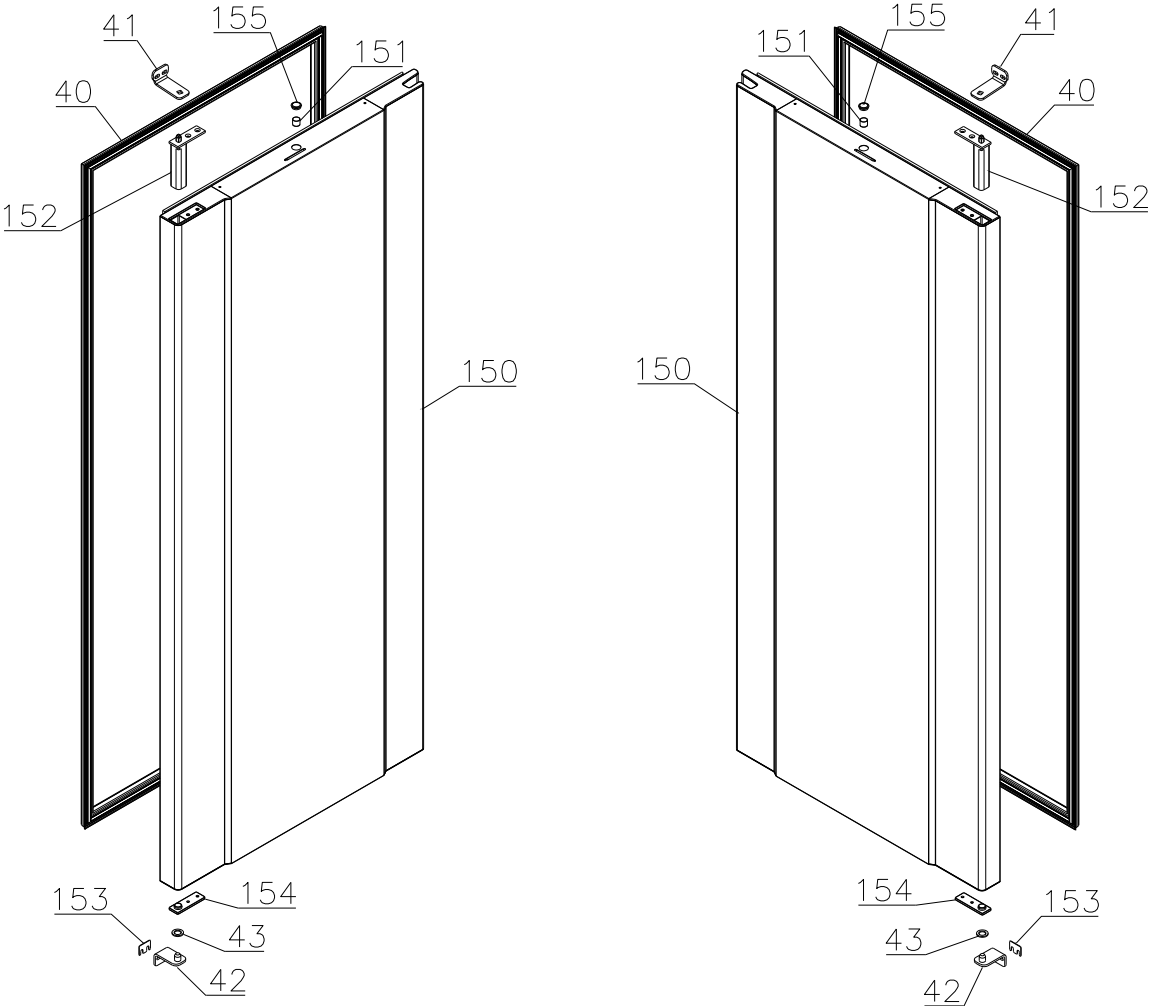


Description :

Spare parts list for BioCompact II
RR 610 G


Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
09.01.09	JP	19.09.25	Jolu	027		

Chart recorder parts

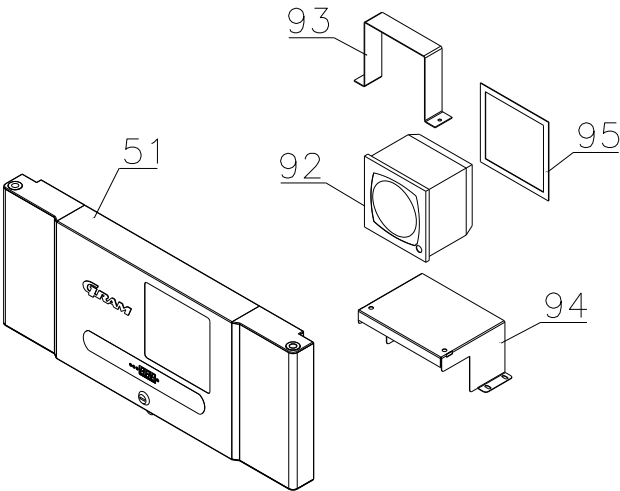

Interior equipment

Shelves:



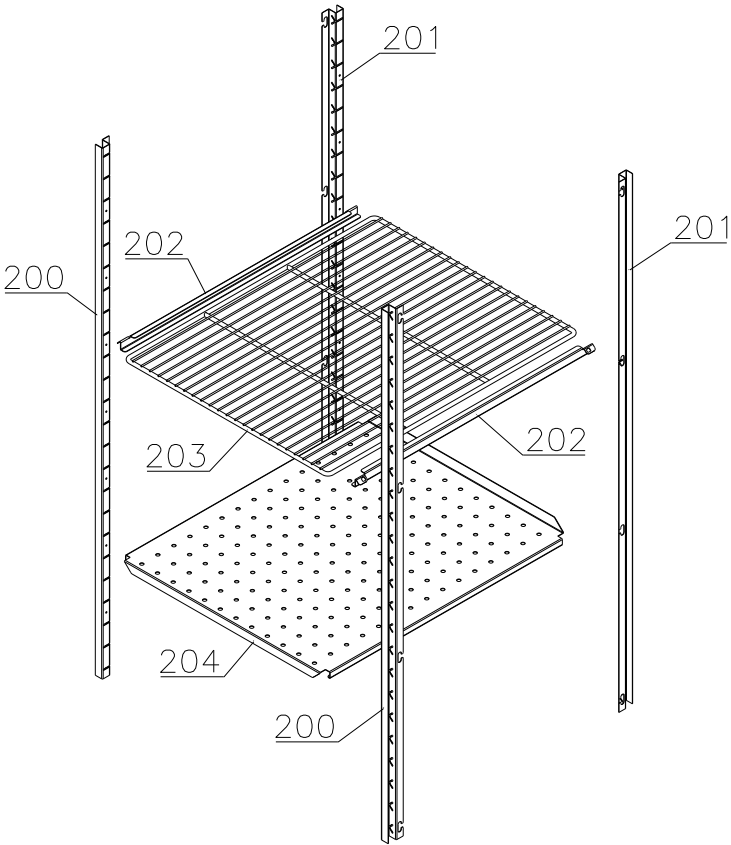


Description :

**Spare parts list for BioCompact II
RR 610 G**

Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
09.01.09	JP	19.09.25	Jolu	027		

Drawers:

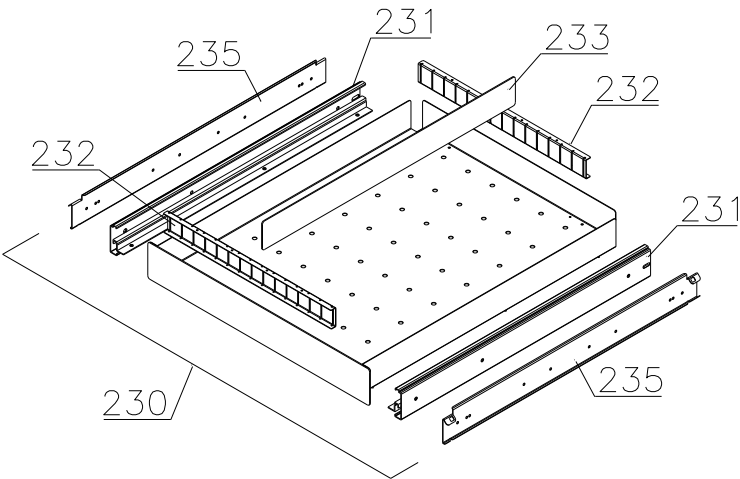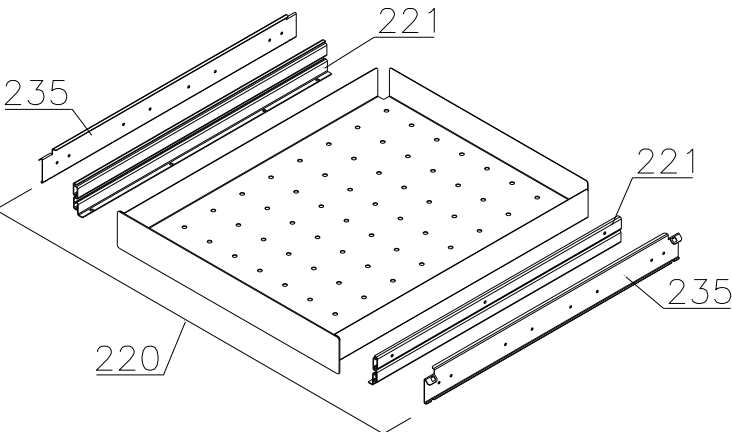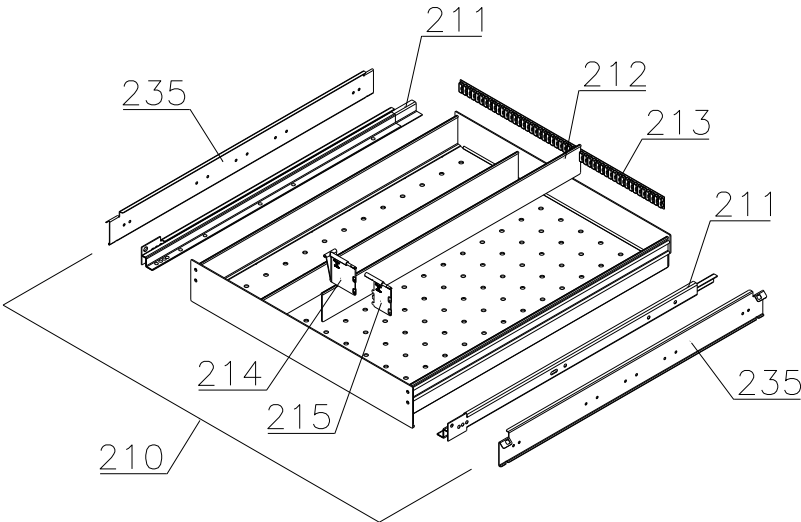

Technical information					765041294	
<u>Description :</u> Spare parts list for BioCompact II RR 610 G						
Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
09.01.09	JP	19.09.25	Jolu	027		

Item no.	Description	Gram part no.	week/year introduced	week/year deleted
----------	-------------	---------------	----------------------	-------------------

Essential spare parts

1	Compressor NLE 11 KK.4	500550024	0508	
2	Starting relay	453050178	0508	
3	Cord relief	452440297	0508	
4	Cover	452440293	0508	
5	Run capacitor 4µF	451723513	0508	
6	Rubber foot	760080519	0508	
7	Screw for compressor	423200210	0508	
8	Fastening clip	760082367	0508	
9	Condenser	612900500	0508	
10	Condenser fan guard compl.	765080189	0508	
11	Filter drier	762900157	0508	
12	Condenser sensor	453531500	0508	
13	Re-evaporation heater	761250778	0508	
14	Float switch	765072049	0508	
15	Evaporator fan complete	761221902	0508	3921
15	Evaporator fan complete	761221904	3921	
16	Evaporator	761220332	0508	
17	Heat exchanger	761221403	0508	
18	Evaporator sensor	453531501	0508	
19	Cabinet sensor	453531501	0508	
20	E-sensor	453531504	0508	
21	Door switch	450400113	0508	
25	Controller display part	612900056	0508	0825
25	Controller display part	612900059	0925	
26	Label with keys	765041752	0508	0825
26	Label with keys	765041889	0925	
27	Band cable	453582046	0508	
27	Band cable for LTP	765072958	3519	
28	Electronic controller K2+	761256120	0508	
28	Electronic controller C01 (EBERHARDT)	761254156	4125	
29	Electrical box complete	761232111	0508	0825
29	Electrical box complete	761233702	0925	
29	Electrical box complete (LTP)	761233700	0925	
30	LTP controller	453580515	3519	
30	EMC/LTP kit - retro fit BioCompact 610	761234073	3519	0925
30	EMC/LTP kit - retro fit BioCompact 610	761234089	1025	
31	LTP sensor	453531511	0718	3519
31	LTP sensor	453531526	3519	
32	PCB protection sheet	765072698	0718	
33	Screening plate	765073271	0718	
34	Dual locking spacer H=7.9 mm	451119525	0718	0825
34	Dual locking spacer H=5.0 mm	451119520	0925	
35	Dual locking spacer H=15.9 mm	451119526	0718	
36	Push spacer H=7.9 mm	451119527	0718	0825
36	Push spacer H=5.0 mm	451119533	0925	

Technical information					765041294	
<u>Description :</u> Spare parts list for BioCompact II RR 610 G						
Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
09.01.09	JP	19.09.25	Jolu	027		

Item no.	Description	Gram part no.	week/year introduced	week/year deleted
37	Solid state relay	450200300	4717	
38	Cable to SSR incl. resistance	765071445	1023	
38	Cable to SSR incl. resistance "LTP"	765071449	1023	
40	Door gasket	761222863	0508	
41	Hinge, top	761220385	0508	4314
41	Hinge, top, self-close right	761222802	4314	
41	Hinge, top, self-close left	761222803	4314	
42	Hinge, bottom	761220387	0508	4314
42	Hinge, bottom, self-close	761222819	4314	
43	Washer	760250524	0508	4314
43	Washer	421708003	4314	

Technical information					765041294	
<u>Description :</u> Spare parts list for BioCompact II RR 610 G						
Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
09.01.09	JP	19.09.25	Jolu	027		

Item no.	Description	Gram part no.	week/year introduced	week/year deleted
----------	-------------	---------------	----------------------	-------------------

Other spare parts

50	Air relief valve	765081206	0508	
51	Front panel	stainless 761231403	0508	0825
51	Front panel	stainless 761234167	0925	
51	Front panel	white 761231404	0508	0825
51	Front panel	white 761234165	0925	
51	Front panel f. chart recorder	stainless 761231451	4410	0825
51	Front panel f. chart recorder	stainless 761234168	0925	
51	Front panel f. chart recorder	white 761231452	4410	0825
51	Front panel f. chart recorder	white 761234166	0925	
52	Hinge for front panel	761220388	0508	
53	Evaporator cover, side/back	761220326	0508	
54	Evaporator cover, front	761220325	0508	
55	Suspension bracket	760250295	0508	
56	Fan holder	761220369	0508	
57	Drip water tray	761221430	0508	
58	Air distribution plate	761220331	0508	
59	Re-evaporation tray	765081149	0508	
60	Water trap	760082891	0508	
61	Drain pipe	765081153	0508	5017
61	Drain pipe	765081253	5017	
62	Heater bracket	761255445	0508	2714
62	Heater bracket	763250002	2714	
63	Insulation slab	761221734	0508	
64	Screw for fan blade	612900100	0508	
65	Float switch bracket	761255446	0508	2714
65	Float switch bracket	763250003	2714	
66	Plastic cap	450450080	0508	
67	Fan guard	765081239	0508	
70	Lead and plug, DK 3-pin	765072568	0508	
70	Lead and plug, Schuko	765072566	0508	
70	Lead and plug, Switzerland	765070163	0508	
70	Lead and plug, GB	765072567	0508	
74	Leg, stainless H=135-200	765080115	0508	
75	Fixed wheel, H=125	480270520	0508	
76	Castor wheel, with brake	480260050	0508	
77	Screw M8 x 12	422715000	0508	
79	Condenser filter	761114416	0508	
80	Attachment spring	762110375	0508	2809
80	Attachment plug	765081246	2809	
81	Front panel close catch	761220375	0508	
82	Front panel close keeper	761253301	0508	
83	Bracket for lock	761220376	0508	
84	Re-evap. heater kit complete	763254026	0508	

<i>Technical information</i>					765041294
Description : Spare parts list for BioCompact II RR 610 G					
Compiled.		Latest revision.			Deleted.
Date	Name	Date	Name	Rev.no	Date
09.01.09	JP	19.09.25	Jolu	027	

Item no.	Description	Gram part no.	week/year introduced	week/year deleted	
85	Door lock	765080092	0508		
86	Key no. 1 (25010)	765080340	0508		
86	Key no. 2 (25020)	765080341	0508		
86	Key no. 3 (25030)	765080342	0508		
86	Key no. 4 (25040)	765080343	0508		
86	Key no. 5 (25050)	765080344	0508		
86	Master key	765080385	0508		
87	Box for alarm connection	761233402	0508		
88	Insulation plug	765080563	0508		
89	Plastic cap	501020023	0508		
90	Liquid container	765080384	0508		
91	Holder for liquid container	761256497	0508		
92	Chart recorder	761256438	4410		
93	Bracket for chart recorder	761256439	4410		
94	Base for chart recorder	761231447	4410		
95	Cover for chart recorder	761256452	4410		
96	Electrics cover	765070143	0508		
97	Set of castor wheels	818899905	0508		
100	Foot pedal complete	819581053	0911		
101	Bracket for pedal	761224010	0911		
102	Pedal arm	761224002	0911		
103	Spindle	761224007	0911		
104	Spring	761224005	0911		
105	Plastic cap	581400021	0911		
106	Bush	761224008	0911		
150	Door complete	stainless	761230917	0508	4314
150	Door complete	white	761230916	0508	4314
150	Door complete, self-close, right	stainless	761232903	4314	4024
150	Door complete, self-close, right	stainless	762234014	4124	
150	Door complete, self-close, left	stainless	761232913	4314	4024
150	Door complete, self-close, left	stainless	762234015	4124	
150	Door complete, self-close, right	white	761232904	4314	4024
150	Door complete, self-close, right	white	762234012	4124	
150	Door complete, self-close, left	white	761232914	4314	4024
150	Door complete, self-close, left	white	762234013	4124	
151	Magnet for door switch	450401002	0508		
152	Self-close device	763251745	4314		
153	Spacer for hinge	761258621	0508		
154	Hinge bearing complete	763251755	4314		
155	Plastic cap, white	760082756	0508		
155	Plastic cap, grey	765081220	0508		

<i>Technical information</i>					765041294
<u>Description :</u> Spare parts list for BioCompact II RR 610 G					
Compiled.		Latest revision.			Deleted.
Date	Name	Date	Name	Rev.no	Date Name
09.01.09	JP	19.09.25	Jolu	027	

Item no.	Description	Gram part no.	week/year introduced	week/year deleted
----------	-------------	---------------	-------------------------	----------------------

Interior equipment:

200	Wall rail, front	761220335	0508	
201	Wall rail, back	761220336	0508	
202	Shelf support rail	761256455	0508	
203	Grey wire shelf	818880008	0508	
204	Perforated shelf, stainless	761230083	0508	
210	Alu drawer, complete	761230362	2210	
211	Set of sliding rails	761234055	2210	
212	Divider rail	761230016	2210	
213	Holder for divider rail	761230063	2210	
214	Plastic divider 1	761230021	2210	
215	Plastic divider 2	761230022	2210	
220	Stainless drawer 30 kg, complete	761231232	0508	
221	Set of sliding rails	761234000	0508	
230	Stainless drawer 40 kg, complete	761231202	0508	
231	Set of sliding rails	761234058	0508	
232	Divider rail	761230305	0508	
233	Divider	761230308	0508	
235	Support for sliding rail	761230201	0508	3310
235	Support for sliding rail, right	761230278	3310	
235	Support for sliding rail, left	761230281	3310	