

*Spare parts -
BioPlus/BioBlood
ER/BR/RF/BF 600/660 W
CH/LH*




Description :

**Spare parts list for BioPlus/BioBlood
ER/BR/RF/BF 600/660 W CH/LH**

Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

ER/BR 600/660 W

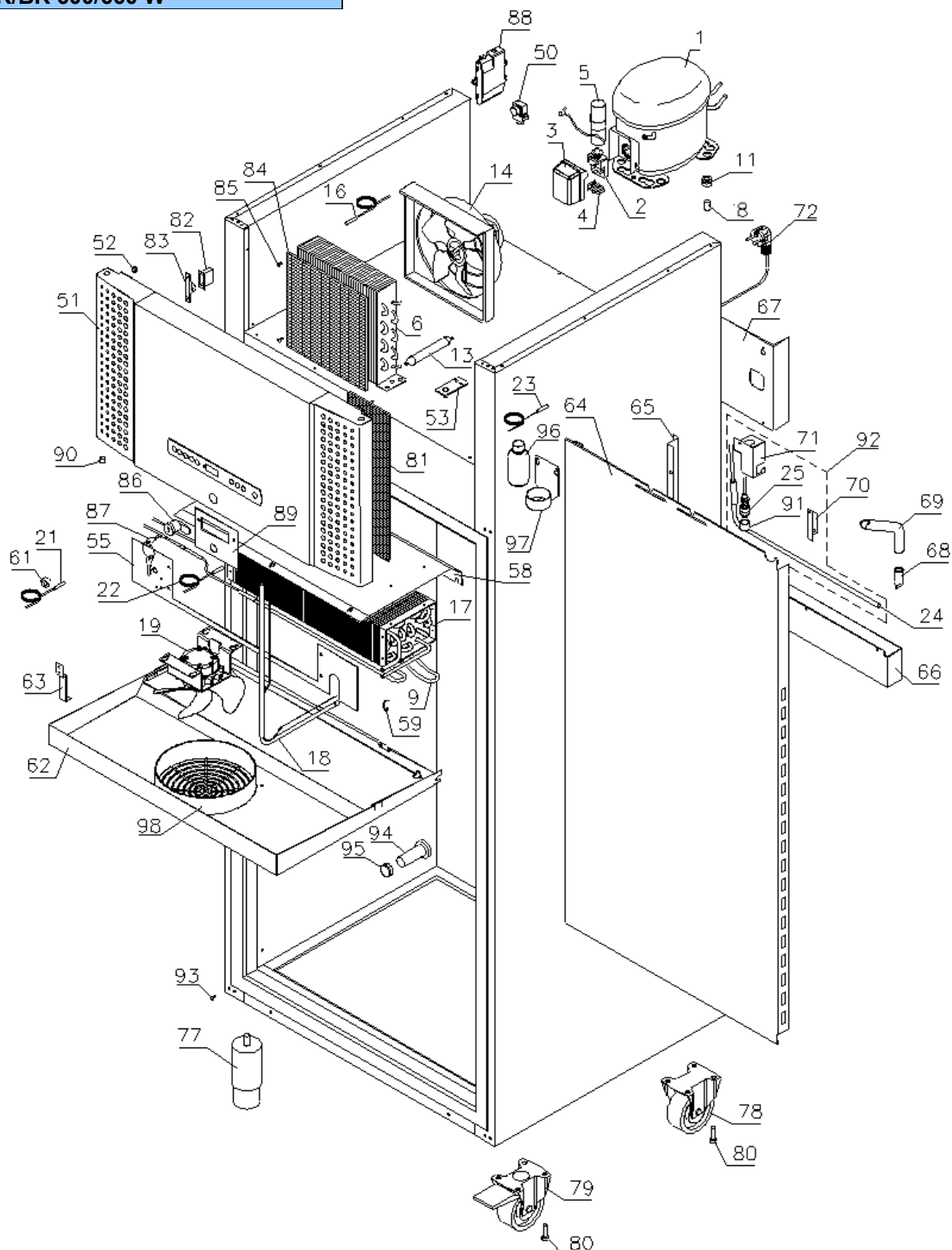


Description :

**Spare parts list for BioPlus/BioBlood
ER/BR/RF/BF 600/660 W CH/LH**

Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

RF/BF 600/660 W– single compressor models

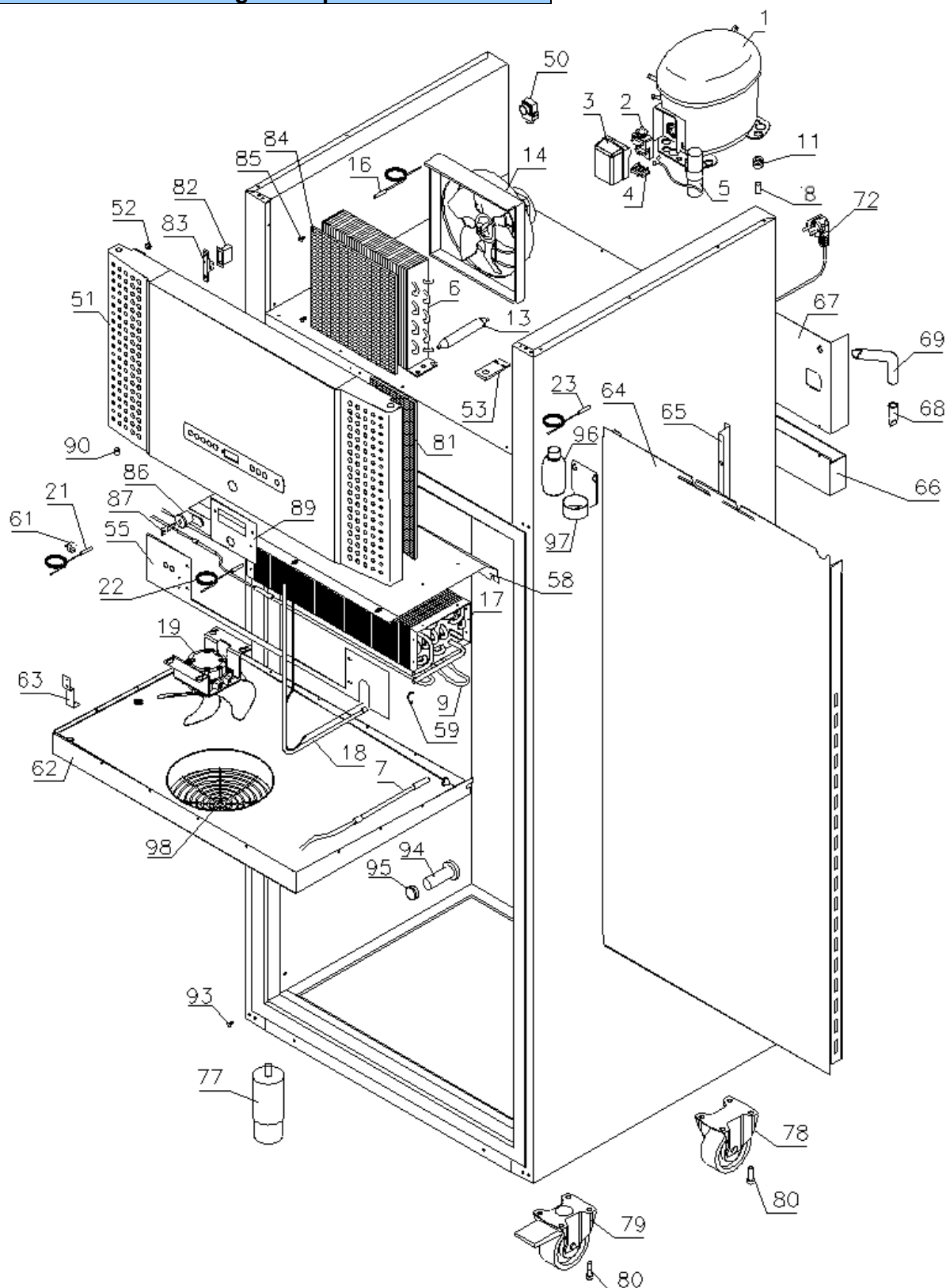


Description :

**Spare parts list for BioPlus/BioBlood
ER/BR/RF/BF 600/660 W CH/LH**

Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

RF/BF 600/660 W – dual compressor models

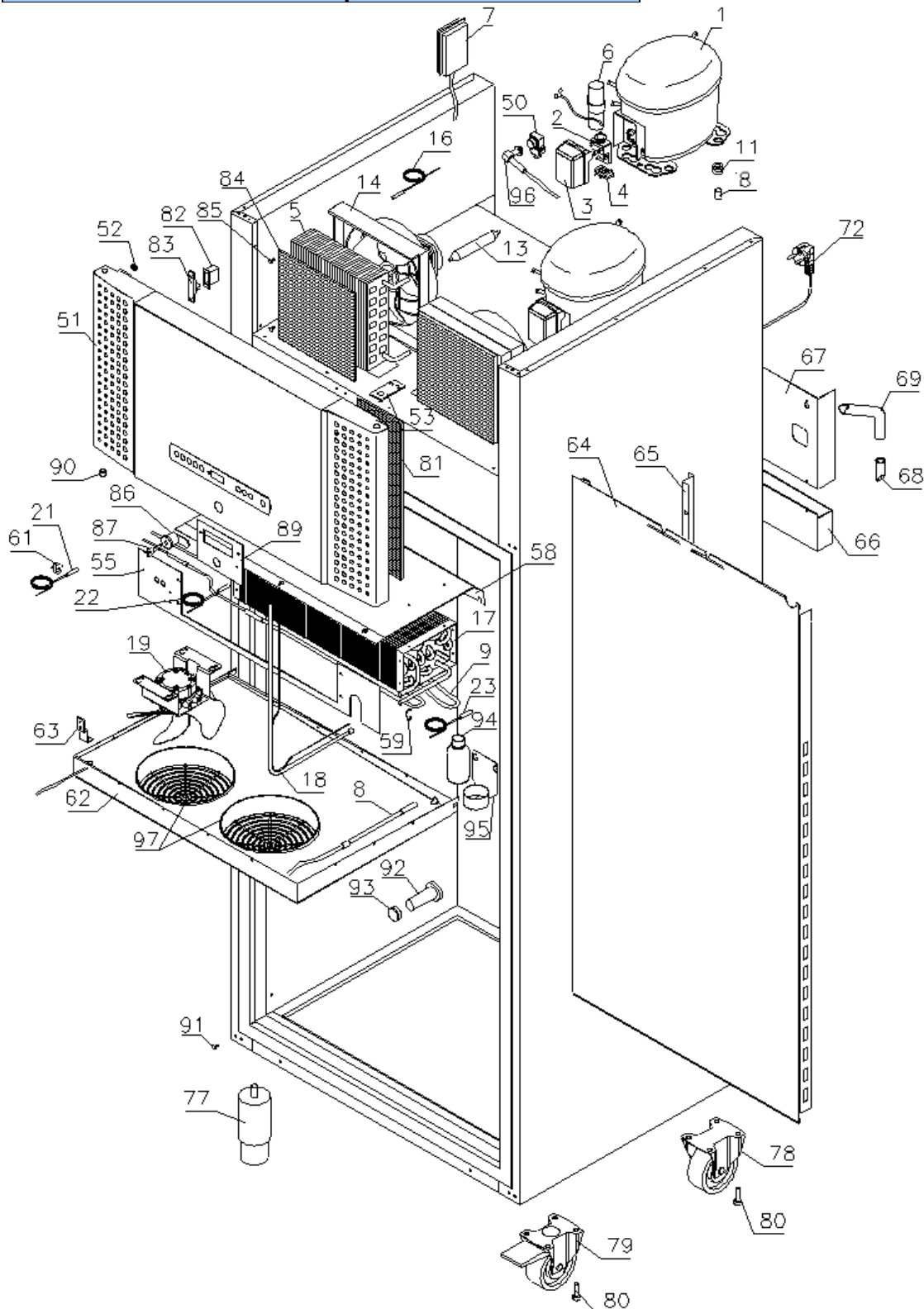


Description :

**Spare parts list for BioPlus/BioBlood
ER/BR/RF/BF 600/660 W CH/LH**


Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

Evaporator cover

1007 – 1009

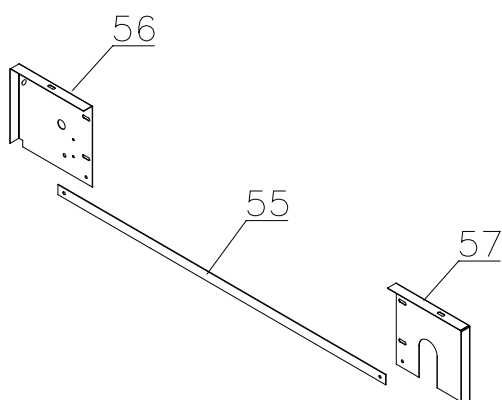


1009 -

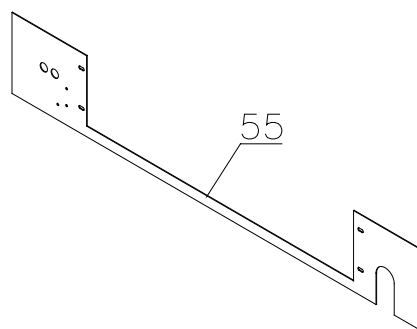


Condenser filter

1007 – 2809

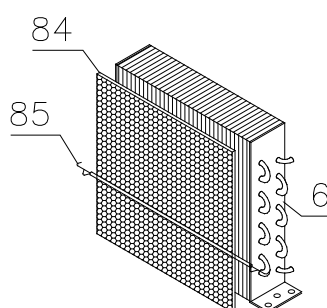

2809 -

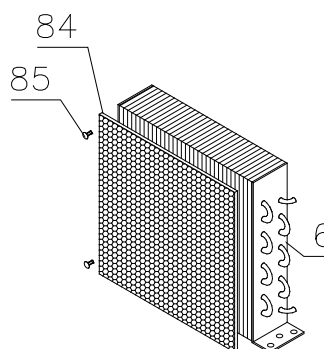

Condenser filter



1007 – 3620

3620 -

Description :

**Spare parts list for BioPlus/BioBlood
ER/BR/RF/BF 600/660 W CH/LH**

Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

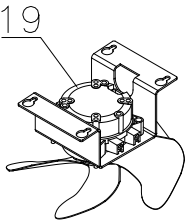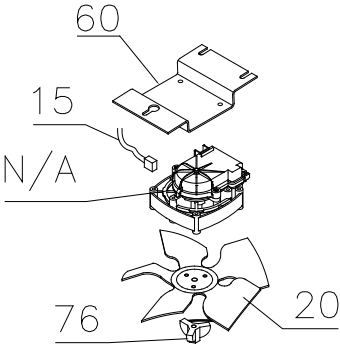


Description :

**Spare parts list for BioPlus/BioBlood
ER/BR/RF/BF 600/660 W CH/LH**


Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

Interior equipment

Shelves:

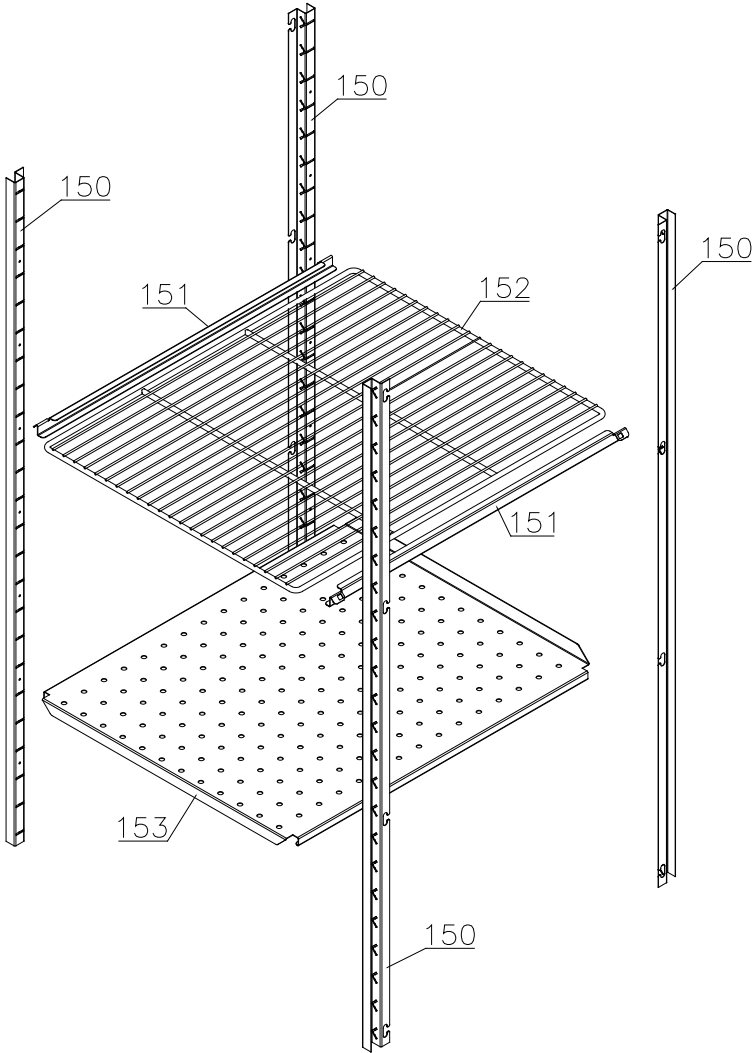

Defrost heater

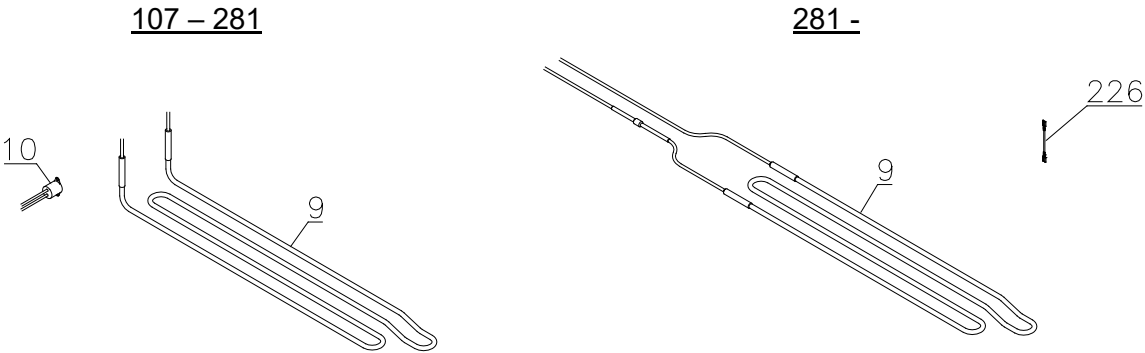


Description :

**Spare parts list for BioPlus/BioBlood
ER/BR/RF/BF 600/660 W CH/LH**


Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

Drawers:


Alu

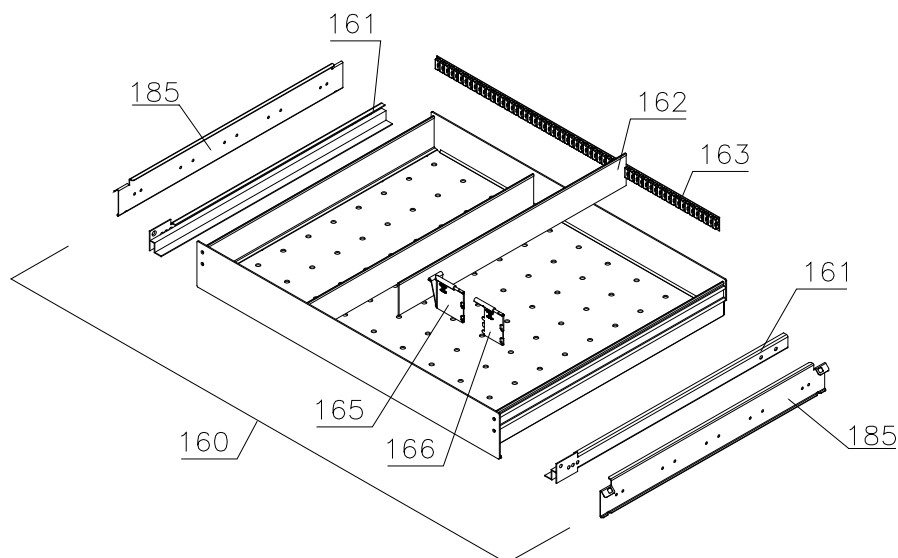

Stainless – 30 kg

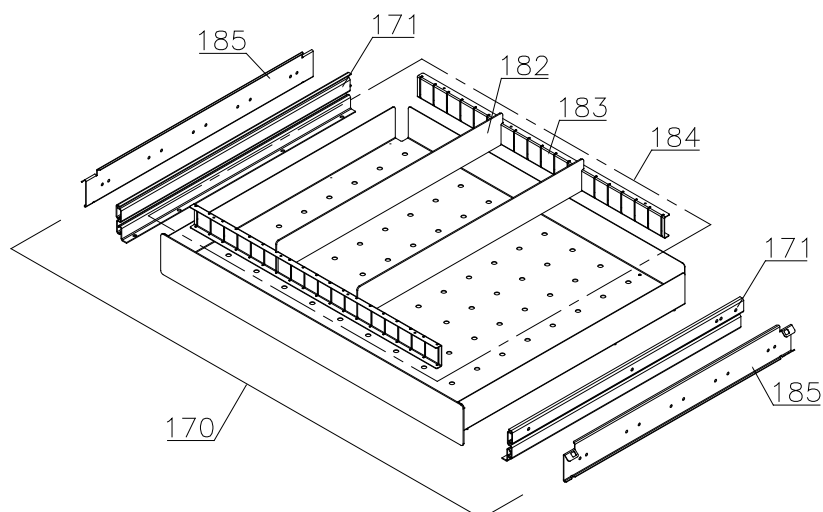

Stainless – 40 kg

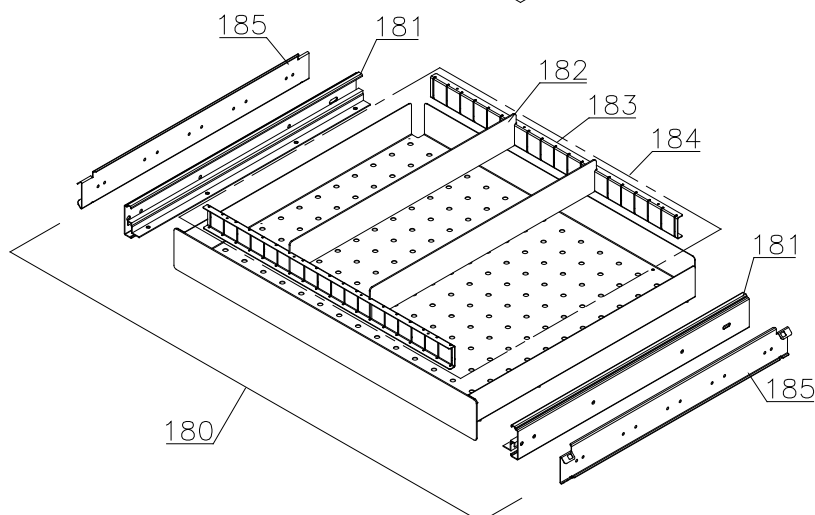


Description :

**Spare parts list for BioPlus/BioBlood
ER/BR/RF/BF 600/660 W CH/LH**


Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

Control panel / electrical parts

1007 - 1709

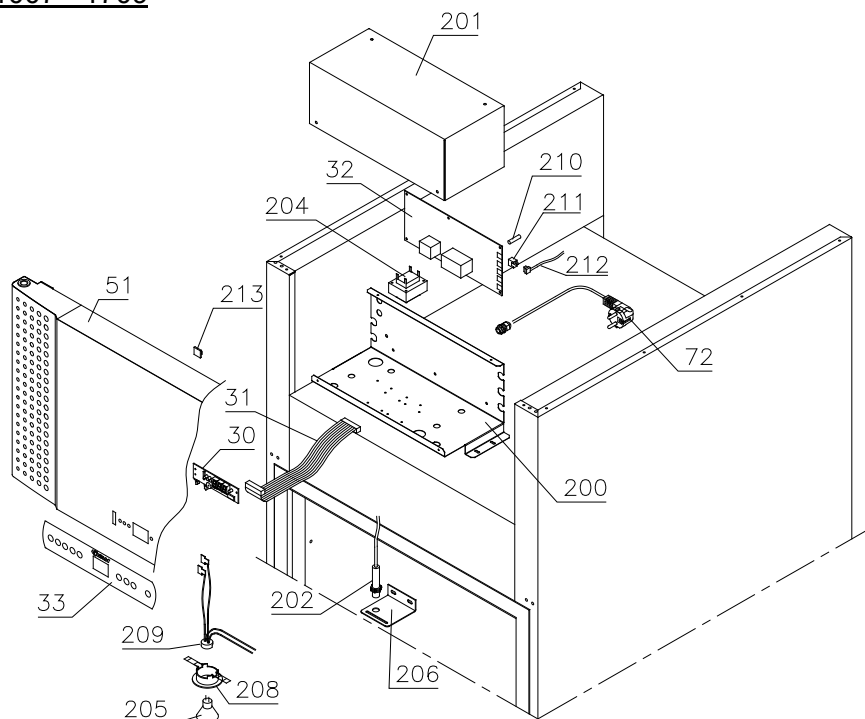

1709 - 4715

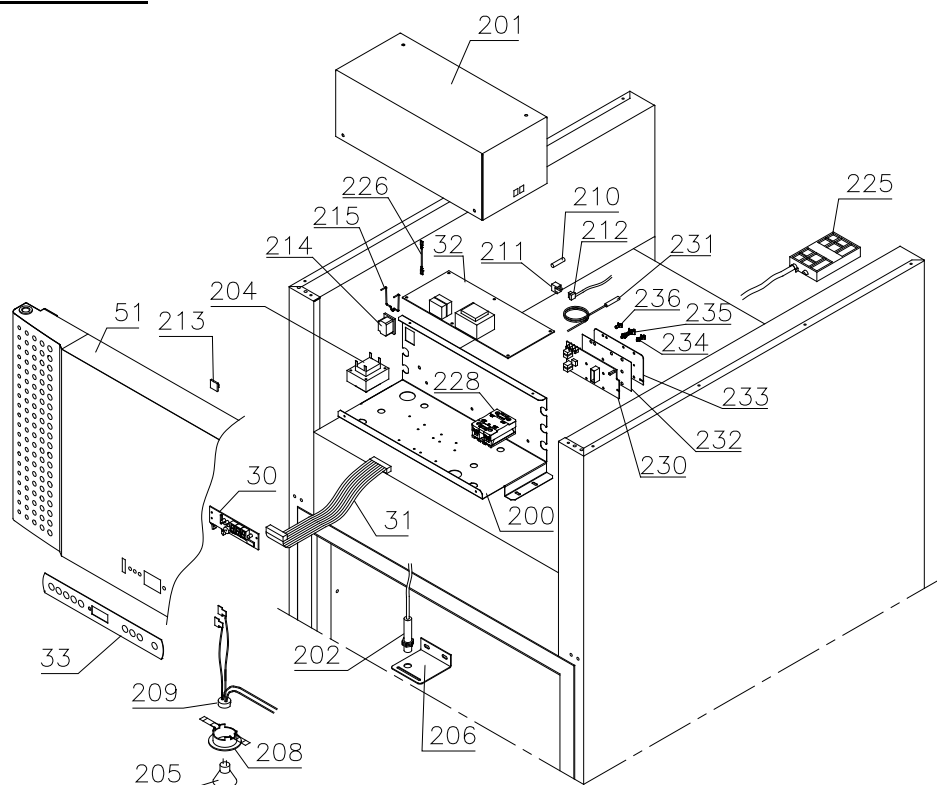


Description :


**Spare parts list for BioPlus/BioBlood
ER/BR/RF/BF 600/660 W CH/LH**

Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

4715 -

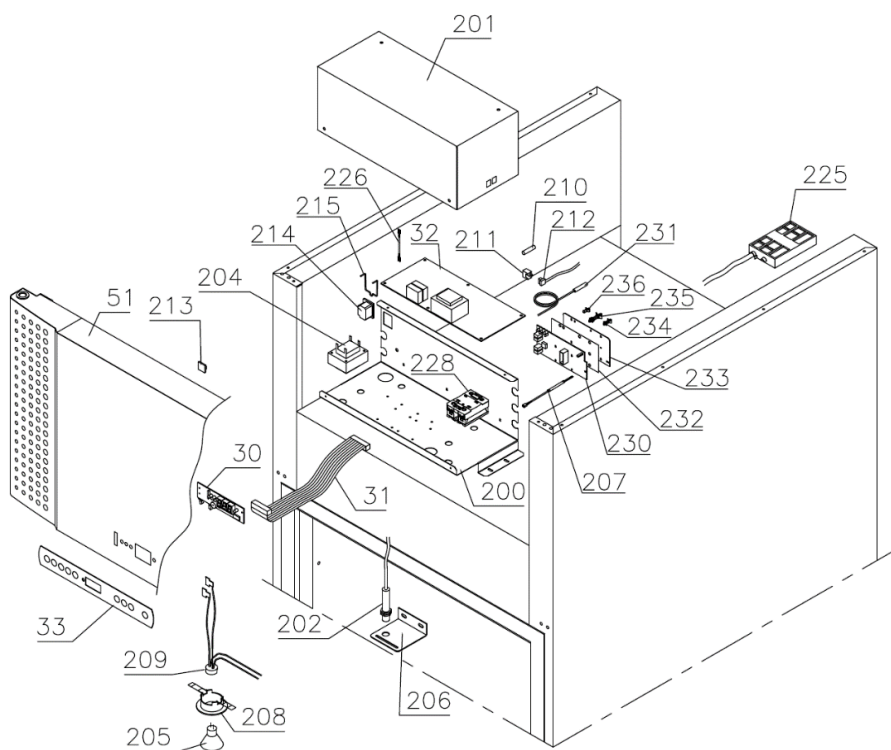

Optional parts

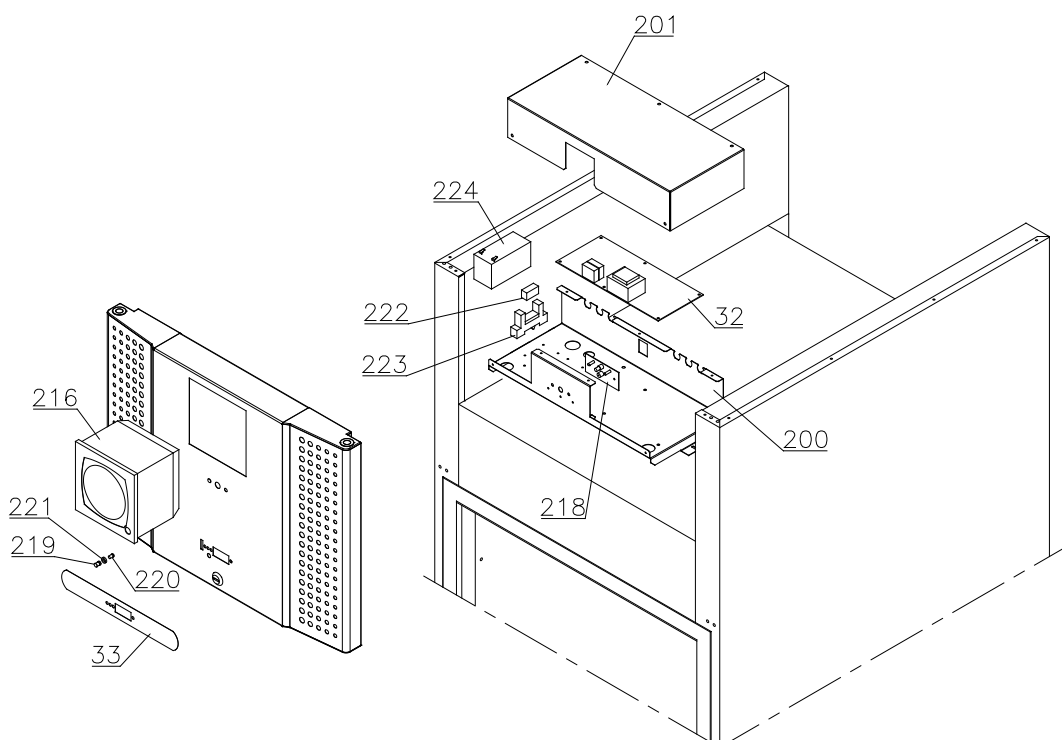


Description :

Spare parts list for BioPlus/BioBlood
ER/BR/RF/BF 600/660 W CH/LH


Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

Door

Lefthand

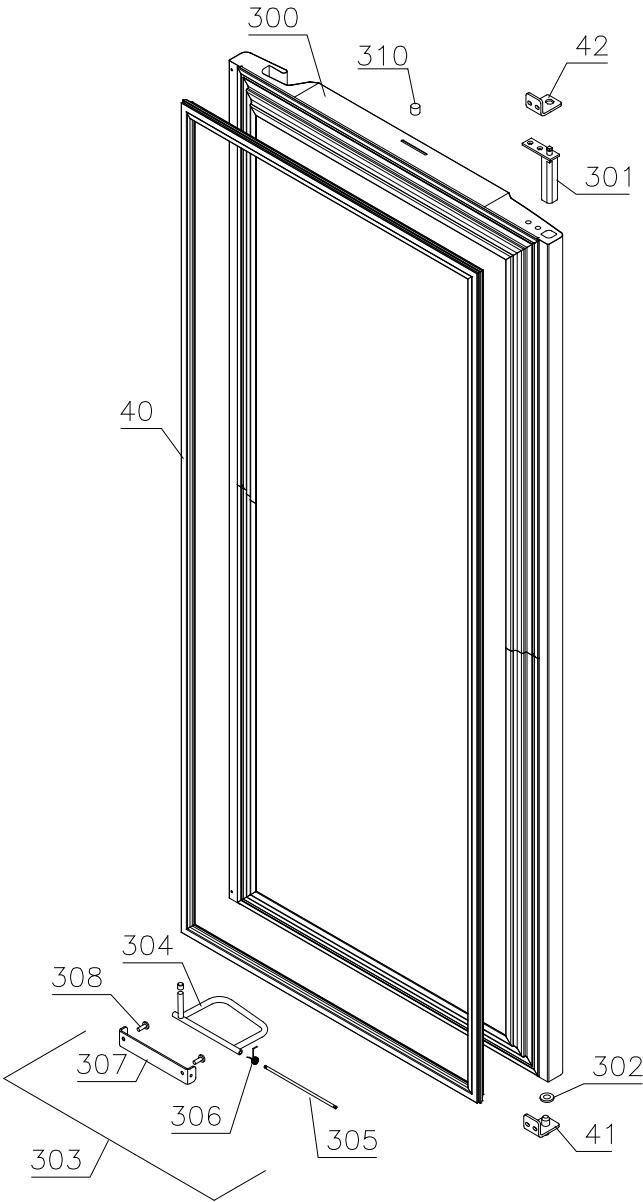

Righthand

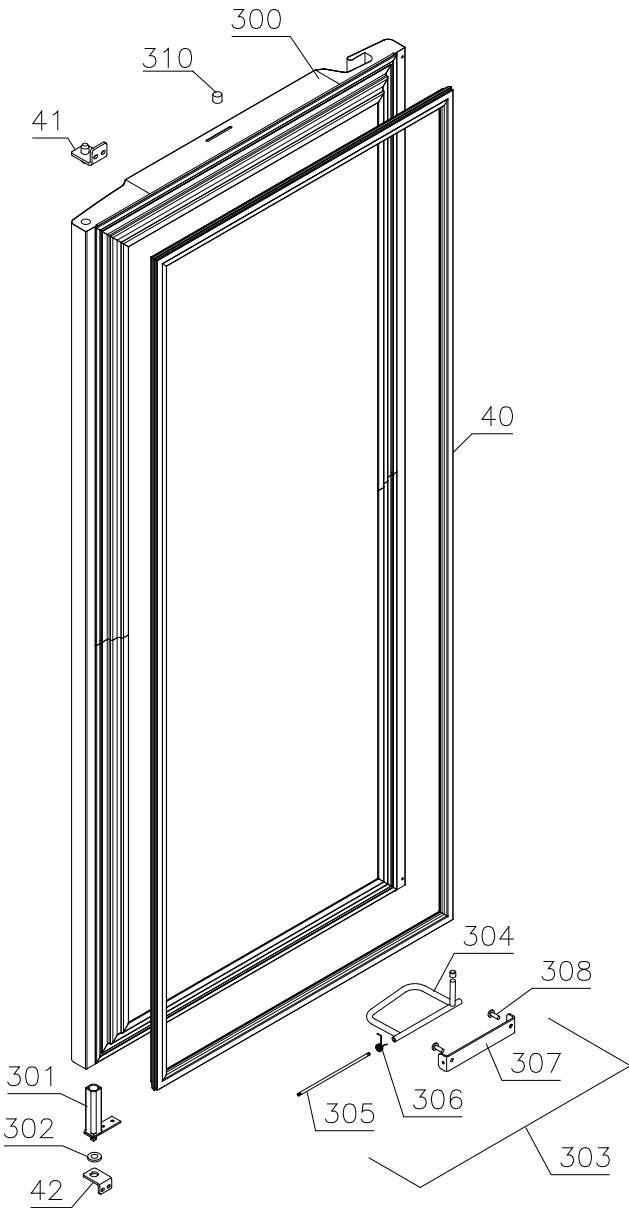


Description :

Spare parts list for BioPlus/BioBlood
ER/BR/RF/BF 600/660 W CH/LH

Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

Glass door

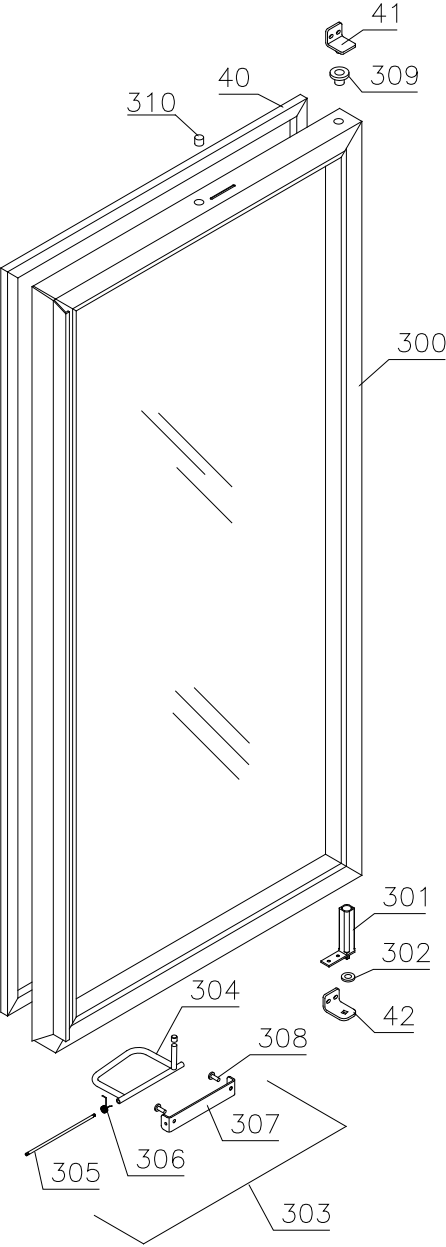


Description :

Spare parts list for BioPlus/BioBlood
ER/BR/RF/BF 600/660 W CH/LH

Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

Internal doors

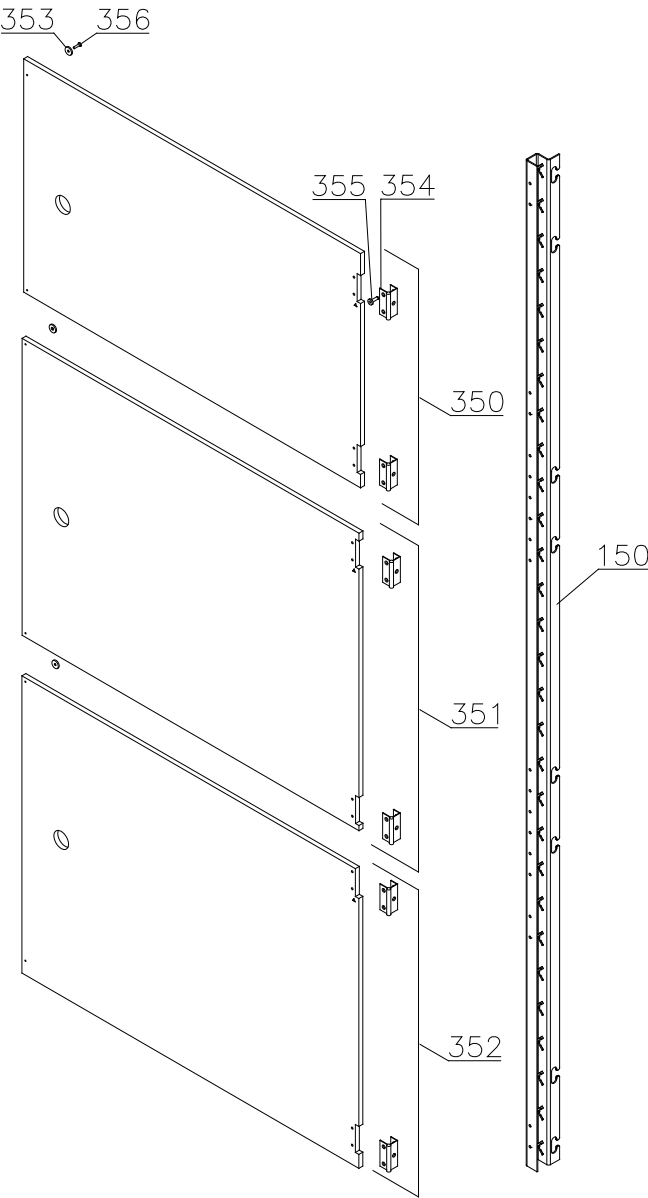

Gram Commercial Technical information					765041221	
<u>Description :</u> Spare parts list for BioPlus/BioBlood ER/BR/RF/BF 600/660 W CH/LH						
Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

Item No.	Description	Gram part No.	week/year Issued	week/year Deleted
----------	-------------	---------------	---------------------	----------------------

Essential spare parts

Refrigeration system/cabinet

ER/BR 600/660 W – 50/60Hz:

1	Compressor FR 7,5 G	500550061	1007
2	Start relay	410589015	1007
3	Cover	452440293	1007
4	Cord relief	452440300	1007
5	Start capacitor	451720018	1007
6	Condenser	765080161	1007

RF/BF 600/660 W – 50Hz:

1	Compressor SC 12 CL	500520255	1007	1613
1	Compressor SC 12 CLX.2	500520257	1613	
2	Start relay	410589038	1007	
3	Cover	452440280	1007	
4	Cord relief	452440292	1007	
5	Start capacitor	451720018	1007	
6	Condenser	765080163	1007	
7	Drip water heater 6,5 W	765071698	1007	

RF/BF 600/660 W – 60Hz:

1	Compressor SC 10 CLX	500520102	1007
2	Start relay	410589021	1007
3	Cover	452440291	1007
4	Cord relief	452440292	1007
5	Start capacitor	451720020	1007
6	Condenser	765080163	1007
7	Drip water heater 6,5 W	765071698	1007

Gram Commercial Technical information					765041221	
<u>Description :</u> Spare parts list for BioPlus/BioBlood ER/BR/RF/BF 600/660 W CH/LH						
Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

Item No.	Description	Gram part No.	week/year Issued	week/year Deleted
----------	-------------	---------------	---------------------	----------------------

Common parts:

8	Compressor screw	760082711	0114	
9	Defrost heating element 350 W	765071819	1007	
10	Overheat safety thermostat	765070828	1007	2811
11	Rubber foot	760080519	1007	
12	Bush	760080521	1007	0313
13	Filter drier	762900156	1007	
14	Condenser fan guard	ER/BR 765080188	1007	
14	Condenser fan guard	RF/BF 765080190	1007	
15	Wire for fan motor	612900400	1007	
16	Condenser sensor	453531500	1007	
17	Evaporator complete	762900090	1007	
17	Evaporator	765080185	1007	
18	Heat exchanger	762900060	1007	
19	Evaporator fan complete	761234098	1007	3921
19	Evaporator fan complete	761234129	3921	
20	Fan blade, ø200	612900201	1007	3620
21	Cabinet sensor	453531500	1007	
22	Defrost sensor	453531501	1007	0712
22	Defrost sensor	453531504	0712	
23	E-sensor	453531504	1007	
24	Re-evap. heating element 60 W	761250778	1007	2714
24	Re-evap. heating element 60 W	765072862	2714	
25	Float switch	765072049	1007	

Gram Commercial Technical information					765041221	
<u>Description :</u> Spare parts list for BioPlus/BioBlood ER/BR/RF/BF 600/660 W CH/LH						
Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

Item No.	Description	Gram part No.	week/year Issued	week/year Deleted
----------	-------------	---------------	---------------------	----------------------

Control panel:

30	Electronic control, display part	612900056	1007	
30	Electronic control, display part for LTP	612900057	3919	
31	Band cable L=1200	453582046	1007	
31	Band cable for LTP	765072969	3919	
32	Controller power part M5	ER 761256135	1007	
32	Controller power part M4	BR 761256133	1007	
32	Controller power part F5	RF/BF 761256134	1007	4414
32	Controller power part F51	RF/BF 761256162	4414	
32	Controller power part M5 batt.alm.	ER 761256146	1007	
32	Controller power part M4 batt.alm.	BR 761256145	1007	
32	Controller power part F5 batt.alm.	RF/BF 761256147	1007	
33	Label with keys	765041752	1007	

Door:

40	Door gasket	600	761259246	1007
40	Door gasket	660	761259247	1007
41	Hinge, upper right/bottom left		761258923	1007
42	Hinge, bottom right/upper left		761258998	1007

Gram Commercial Technical information					765041221	
<u>Description :</u> Spare parts list for BioPlus/BioBlood ER/BR/RF/BF 600/660 W CH/LH						
Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

Item No.	Description	Gram part No.	week/year Issued	week/year Deleted
----------	-------------	---------------	---------------------	----------------------

Other spare parts

Cabinet

50	Air relief valve		765081206	1007	
51	Front panel – solid door, stainless		761231464	1007	
51	Front panel – solid door, white		761231465	1007	
51	Front panel – glass door, stainless		761231462	1007	
51	Front panel – glass door, white		761234025	1007	
52	Front panel plug		760250272	1007	
53	Hinge for front panel		761258926	1007	
55	Cover rail for evaporator		761254449	1007	1009
55	Cover for evaporator (one piece)		761254469	1009	
56	Cover plate, left		761255454	1007	1009
57	Cover plate, right		761255453	1007	1009
58	Suspension plate for evaporator		761255455	1007	
59	Heating element bracket		761254427	1007	
60	Fan holder		761258499	1007	3620
61	Cable clip		452470001	1007	
62	Drip water tray		761256013	1007	5108
62	Drip water tray		761256032	5108	
62	Drip water tray, insulated Single	RF/BF	761256048	0700	
62	Drip water tray, insulated Dual	RF/BF	761256044	0700	
63	Holder for drip water tray		761253441	1007	
63	Holder for drip water tray	RF/BR	761253442	0710	
64	Air distribution plate	600	761254480	1007	
64	Air distribution plate	660	761254481	1007	
65	Rail for air distribution plate		760243634	1007	
66	Re-evaporation tray		765081149	1007	
67	Cover plate for re-evap. tray		761255383	1007	
68	Water trap		760082891	1007	
69	Drain pipe		765081253	1007	
70	Heating element bracket		761255445	1007	2714
70	Heating element bracket		763250002	2714	
71	Float switch bracket		761255446	1007	2714
71	Float switch bracket		763250003	2714	
72	Lead and plug, DK 3-pin		765072555	1007	
72	Lead and plug, Schuko		765072554	1007	
72	Lead and plug, GB		765072553	1007	

Gram Commercial Technical information					765041221	
<u>Description :</u> Spare parts list for BioPlus/BioBlood ER/BR/RF/BF 600/660 W CH/LH						
Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

Item No.	Description	Gram part No.	week/year Issued	week/year Deleted
76	Screw for fan blade	612900100	1007	3620
77	Leg, stainless H=135-200	765080115	1007	
78	Fixed wheel, H=125	480270520	1007	
79	Castor wheel, with brake	480260050	1007	
80	Screw M8 x 12	420050100	1007	
81	Front panel filter	761258491	1007	
82	Front panel close latch	761253307	1007	
83	Front panel close keeper	761253301	1007	
84	Condenser filter	761114415	1007	
85	Attachment spring	761114417	1007	2809
85	Attachment plug	765081233	2809	
86	Door lock – solid door	765080092	1007	
86	Door lock – glass door	765080098	1007	
87	Key no. 1 (25010)	765080340	1007	
87	Key no. 2 (25020)	765080341	1007	
87	Key no. 3 (25030)	765080342	1007	
87	Key no. 4 (25040)	765080343	1007	
87	Key no. 5 (25050)	765080344	1007	
87	Master key	765080385	1007	
88	Junction box f. re-evap. heater	765070143	1007	
89	Bracket for lock	761258747	1007	
90	Bush ø12	761254345	1007	
91	Plastic cap	450450080	1007	
92	Re-evaporation heater kit complete	761255438	1007	2714
92	Re-evaporation heater kit complete	763254026	2714	
93	Cap for M5 hole grey	765080501	1007	
93	Cap for M5 hole white	765080504	1007	
94	Insulation plug	765080563	2309	
95	Plastic cap	501020023	2309	
96	Liquid container	765080384	1007	
97	Holder for liquid container	761256497	1007	
98	Fan guard	765081239	0509	

Gram Commercial Technical information					765041221	
<u>Description :</u> Spare parts list for BioPlus/BioBlood ER/BR/RF/BF 600/660 W CH/LH						
Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

Item No.	Description	Gram part No.	week/year Issued	week/year Deleted
----------	-------------	---------------	---------------------	----------------------

Interior equipment:

150	Wall rail	600	761256462	1007
150	Wall rail - for internal doors	600	761230295	2010
150	Wall rail	660	761256454	1007
150	Wall rail - for internal doors	660	761230296	2010
151	Shelf support rail (single piece)		760248857	1007
151	Shelf support rail (set of 2)		818790030	1007
152	Stainless wire shelf		818799967	1007
152	Stainless wire shelf / internal doors		765080403	3210
152	Grey wire shelf		818799968	1007
153	Perforated shelf, stainless		761230058	1007
160	Alu drawer, complete		761230358	1007
161	Set of sliding rails		761234056	1007
162	Divider rail		761230053	1007
163	Holder for divider rail		761230366	1007
165	Plastic divider 1		761230021	1007
166	Plastic divider 2		761230022	1007
170	Stainless drawer, complete 30 kg		761231233	1007
171	Set of sliding rails		761234007	1007
180	BioBlood drawer, complete 40 kg		761231200	1008
181	Set of sliding rails		761234057	1008
182	Divider, stainless		761230312	2409
183	Holder for divider, stainless		761230311	2409
184	Divider set, complete		761231449	2409
185	Support for sliding rail		761230218	1008

Gram Commercial Technical information					765041221	
<u>Description :</u> Spare parts list for BioPlus/BioBlood ER/BR/RF/BF 600/660 W CH/LH						
Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

Item No.	Description	Gram part No.	week/year Issued	week/year Deleted
----------	-------------	---------------	---------------------	----------------------

Control panel/electrical parts:

200	Junction box	761230090	1007
200	Junction box - battery alarm	761230089	1007
201	Cover for junction box	761255432	1007
201	Cover for junction box – battery alarm	761230175	1007
202	Door switch	450400113	1007
203	Distance tube L=8.3 mm	765080021	1007
204	Trafo for lights	760073650	1007
205	Halogen light 10W	450742002	1007
206	Bracket for door switch and lock	761259005	1007
207	Cable to SSR incl. resistance	765071445	1023
207	Cable to SSR incl. resistance “LTP”	765071449	1023
208	Mounting socket for halogen light	761251323	1007
209	Wiring for halogen light	765071603	1007
210	Spacer for PCB	451119515	1007
211	Sensor plug holder	452260601	1007
212	Sensor plug	452260600	1007
213	Cable clip	424350001	1007
214	Appliance inlet	451022002	1709
215	Brace for mains cable	761230215	1709
216	Chart recorder	761256438	1007
218	Battery alarm PCB	453580116	1007
219	Clear glass for LED	765080380	1007
220	Light stick for LED	765080381	1007
221	Spacer for LED	765080382	1007
222	Relay	450200082	1007
223	Relay socket	450200097	1007
224	Battery	130080060	1007
225	Box for alarm connection	761230356	1210
226	Wire for klaxon	765072664	2811
227	Datalogger connector	453570037	4311
230	LTP controller	453580515	3919
231	LTP sensor	453531530	3919
231	LTP sensor	453531533	3919
232	PCB protection sheet	765072698	0718
233	Screening plate	765073271	0718
234	Dual locking spacer H=7.9 mm	451119525	0718
235	Dual locking spacer H=15.9 mm	451119526	0718
236	Push spacer H=7.9 mm	451119527	0718

Gram Commercial Technical information					765041221	
<u>Description :</u> Spare parts list for BioPlus/BioBlood ER/BR/RF/BF 600/660 W CH/LH						
Compiled.		Latest revision.			Discontinued	
Date	Name	Date	Name	Rev.no	Date	Name
29.05.07	JP	20.09.24	JOLU	037		

Item No.	Description	Gram part No.	week/year Issued	week/year Deleted
----------	-------------	---------------	---------------------	----------------------

Door :

300	Door, stainless	600	761251303	1007
300	Door, stainless	660	761251308	1007
300	Door, white	600	761251305	1007
300	Door, white	660	761251304	1007
300	Glass door	600	761259169	1007
300	Glass door	660	761259167	1007
301	Self-close device		761259909	1007
302	Washer, righthand		421708020	1007
302	Washer, lefthand		760250524	1007
303	Foot pedal complete, lefthand		761258167	1007
303	Foot pedal complete, righthand		761258166	1007
304	Pedal arm, lefthand		761258482	1007
304	Pedal arm, righthand		761258407	1007
305	Pedal spindle		761258406	1007
306	Spring for pedal, lefthand		760251364	1007
306	Spring for pedal, righthand		760251363	1007
307	Bracket for pedal		761258404	1007
308	Screw M5x12		420178203	1007
309	Hinge bush		760242041	1007
310	Magnet for door switch		450401002	1007

Internal doors:

350	Upper door, complete		761231889	2010
351	Intermediate door, complete	600	761231890	2010
351	Intermediate door, complete	660	761231891	2010
352	Lower door, complete		761231892	2010
353	Magnet		450401003	2010
354	Hinge		761230343	2010
355	Screw for hinge		420168174	2010
356	Screw for magnet		420168175	2010